


Anti Rat Dynein Polyclonal Antibody

Package Size	25 μ g (100 μ L / vial)
Format	Rabbit polyclonal antibody 0.25mg/mL
Buffer	PBS [containing 2% Block Ace as a stabilizer, 0.1% Proclin as a bacteriostat]
Storage	Store below -20°C Once thawed, store at 4°C. Repeated freeze-thaw cycles should be avoided.
Purification method	This antibody was purified from rabbit serum immunized with synthesized partial peptide of rat Dynein by peptide affinity chromatography.
Working dilution	For western blotting ; 0.5~2 μ g/mL

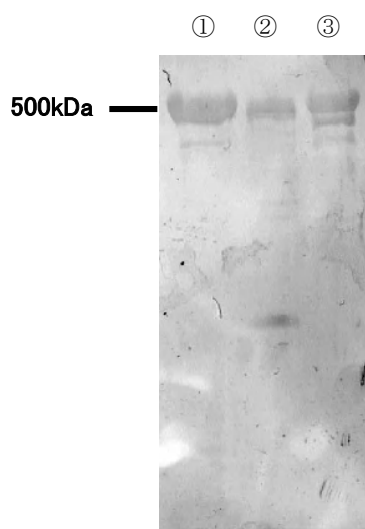

Western Blotting

Sample :

- ① Dynein from pig brain cytoplasmic
- ② 14S Dynein from *Tetrahymena* Axonemes
- ③ 22S Dynein from *Tetrahymena* Axonemes

Preparation of antibodies and instruction :

Toyoshima YY. and Kobayashi T.

Department of Life Sciences,

Graduate School of Art and Science, The University of Tokyo

Anti Rat Dynein Polyclonal Antibody**【Reference】**

1. Gibbons I.R.:Proc. Natl. Acad. Sci. USA 1963,50,1002-1010
2. Hirokawa E, et al.: Proc Natl Acad Sci USA 2000 Mar 14;97(6):2533-7
3. Shirogushi K,et al.: Cell Motil Cytoskeleton 2002 Aug;49(4):189-99
4. Toyoshima YY.: J Cell Biol 1987 Aug; 105(2):897-901
5. Toyoshima YY.: J Cell Biol 1987 Aug; 105(2):887-895
6. Vale RD, et al.: Cell 1988 Feb 12;52(3):459-69
7. Shimizu T,et al.: Biochemistry 1995 Feb 7;34(5):1575-82
8. Yu H, et al.: J Biol Chem 1992 Oct 5;267(28):20457-64
9. Lacey ML,et al.: J Biol Chem 1992 Mar 5;267(7):4793-8

Supplier**SCETI**
SCETI K.K.

3-6-7 Kasumigaseki, Chiyoda-ku, Tokyo 100-0013, JAPAN
Tel +81(3) 5510-2347 Fax +81(3) 5510-0133
URL: <http://www.sceti.jp/export/> e-mail: exp-pet@sceti.co.jp