



Peninsula Laboratories, LLC

A Member of the Bachem Group

305 Old County Road, San Carlos, CA 94070

Tel: (800) 922-1516 • (650) 592-5392

Fax: (650) 595-4071

www.bachem.com

Royalmount Aqueous Mounting Medium

Royalmount is an aqueous mounting medium for cover-slipping tissue sections stained with alcohol soluble chromogens like fast red or AEC (3-amino-9-ethylcarbazole).

Product Number:	T-3402
Quantity:	60ml
	Supplied as liquid solution ready to use
Stability:	Original vial: 1 year at room temperature
Applications:	Before using Royalmount remove the red cap of the bottle and cut off the tip of the nozzle with a sharp blade or a scissors. Make sure that the opening is not too wide. Do not shake the bottle to avoid air bubbles.

Typical Protocol:

- Remove excess liquid from the slide. Tissue must still be humid.
- Apply 1-3 drops of Royalmount to the tissue section.
- Place slide horizontally on a heating plate at 60°C for at least 30 minutes.
- Remove slide from the heating plate and allow it to reach room temperature.
- To remove Royalmount, soak slide overnight in water.
- If required, post-mounting a coverslip on the hardened Royalmount is possible. Apply a few drops of a commercial mounting medium for coverslips on the slide and mount the coverslip.

Royalmount is also available in 250ml size (product T-3403)

For in vitro research only.