



Peninsula Laboratories, LLC

A Member of the Bachem Group

305 Old County Road, San Carlos, CA 94070

Tel: (800) 922-1516 • (650) 592-5392

Fax: (650) 595-4071

www.bachem.com

Cat. No. T-4577 Normal Rabbit Serum (Pretitered)

This serum has been designed to be used specifically in conjunction with pretitered Goat Anti-Rabbit IgG (GARGG) for precipitation of "bound" counts in the titration of antisera or in RIAs.

Reconstitution Information:

Rehydrate the lyophilized powder with 50ml of buffer, this should be the same buffer used in the assays.

Recommended Use:

After the addition of 100ul of GARGG to each tube, add 100ul of the rehydrated NRS to each tube and vortex the contents. Allow the precipitates to form for 2 hours at room temperature. Centrifuge the tubes at 1700xg for 20 minutes and then carefully aspirate off the supernatant. Count the pellet in a scintillation well gamma counter.