



Peninsula Laboratories, LLC

A Member of the Bachem Group

305 Old County Road, San Carlos, CA 94070

Tel: (800) 922-1516 • (650) 592-5392

Fax: (650) 595-4071

www.bachem.com

Cat. No. T-5051

Goat anti-Guinea Pig IgG Serum (Pretitered)

This antiserum has been titered against normal Guinea Pig serum and is designed to be used specifically in conjunction with pretitered normal rabbit serum (catalog #NGS-500) for precipitation of "bound" counts in the titration of antisera or in RIAs.

Reconstitution Information:

Rehydrate the lyophilized powder with 50ml of buffer, this should be the same buffer used in the assays.

Recommended Use:

Add 100ul of the rehydrated GAGGG to each tube, then add 100ul of pretitered normal Guinea Pig serum to each tube and vortex mix the contents. Allow the precipitates to form for 2 hours at room temperature or 6 hours at 4°C. Centrifuge the tube at 1700xg for 20 minutes to collect the precipitates and carefully aspirate off the supernatant. Count the pellet in a scintillation well gamma counter.