



## **Peninsula Laboratories, LLC**

**A Member of the Bachem Group**

305 Old County Road, San Carlos, CA 94070

Tel: (800) 922-1516 • (650) 592-5392

Fax: (650) 595-4071

[www.bachem.com](http://www.bachem.com)

**Cat. No. T-5052**

### **Normal Guinea Pig Serum (Pretitered)**

This serum has been designed to be used specifically in conjunction with pretitered Goat anti-Guinea Pig IgG (GAGGG) for precipitation of "bound" counts in the titration of antisera or in RIA's.

#### **Reconstitution Information:**

Rehydrate the lyophilized powder with 50ml of buffer, this should be the same buffer used in the assays.

#### **Recommended Use:**

After the addition of 100 $\mu$ l of GAGGG to each tube, add 100 $\mu$ l of the rehydrated NGS to each tube and vortex the contents. Allow the precipitates to form for 2 hours at room temperature. Centrifuge the tubes at 1700xg for 20 minutes and then carefully aspirate off the supernatant. Count the pellet in a scintillation well gamma counter.