



## **Peninsula Laboratories, LLC**

**A Member of the Bachem Group**

305 Old County Road, San Carlos, CA 94070

Tel: (800) 922-1516 • (650) 592-5392

Fax: (650) 595-4071

[www.bachem.com](http://www.bachem.com)

**Cat. No. Y-1050**

**RIA Buffer**

### **Reconstitution Information:**

Each vial of the RIA buffer comes with 50 ml of concentrated RIA buffer. Dilute the buffer to a total volume of 200 ml (4x concentration) to be used for 125 tests in our RIA system.

### **Recommended Use:**

This buffer can be used to dilute the components in our RIA kit. For use with 500 test pre-titrated antiserum, 4 x 125 test vials will be provided. After 4-fold dilution, the final working concentration contains 0.1% Triton X-100, 0.1% BSA, 0.01% sodium azide ( $\text{NaN}_3$ ), 0.05M NaCl, and 0.1M sodium phosphate buffer (pH 7.4).