

sHLA-G ELISA

Product Data Sheet

Cat. No.: RD194070100R

For Research Use Only

CONTENTS

1.	INTENDED USE	3
2.	STORAGE, EXPIRATION	3
3.	INTRODUCTION	4
4.	TEST PRINCIPLE	4
5.	PRECAUTIONS	5
6.	TECHNICAL HINTS	5
7.	REAGENT SUPPLIED	6
8.	MATERIAL REQUIRED BUT NOT SUPPLIED	6
9.	PREPARATION OF REAGENTS	7
10.	PREPARATION OF SAMPLES	8
11.	ASSAY PROCEDURE	9
12.	CALCULATIONS	11
13.	PERFORMANCE CHARACTERISTICS	12
14.	DEFINITION OF THE CALIBRATOR	14
15.	METHOD COMPARISON	14
16.	TROUBLESHOOTING AND FAQs	15
17.	REFERENCES	16
18.	EXPLANATION OF SYMBOLS	18

**»» This kit is manufactured in cooperation between:
BioVendor – Laboratorní medicína a.s. and EXBIO Praha a.s.**

»» Use only the current version of Product Data Sheet enclosed with the kit!

1. INTENDED USE

The RD194070100R sHLA-G ELISA is a sandwich enzyme immunoassay for the quantitative measurement of soluble forms of human leukocyte antigen-G (sHLA-G).

»» Features

- **It is intended for research use only**
- The total assay time is about 20 hours
- The kit measures shedded HLA-G1 and HLA-G5 in serum, plasma, amniotic fluid or cell culture supernatant
- Calibrator is human native protein
- Assay format is 96 wells
- Components of the kit are provided ready to use, concentrated or lyophilized

2. STORAGE, EXPIRATION

Store the complete kit at 2-8°C. Under these conditions, the kit is stable until the expiration date (see label on the box).

For stability of opened reagents see Chapter 9.

3. INTRODUCTION

Human leukocyte antigen-G (HLA-G) differs from the other MHC class I genes by its low polymorphism and alternative splicing that generates seven HLA-G proteins, whose tissue-distribution is restricted to normal fetal and adult tissues that display a tolerogenic function toward both innate and acquired immune cells. Soluble HLA-G is an immunosuppressive molecule inducing apoptosis of activated CD8(+) T cells and down-modulating CD4(+) T cell proliferation.

Recently, using specific ELISA to analyse the presence of sHLA-G molecules in culture supernatants of early embryos obtained by in vitro fertilization (IVF) before transfer, several reports demonstrated that positive embryo implantations occurred with embryos secreting sHLA-G molecules. These breakthrough results indicate that sHLA-G ELISA can be a useful biochemical assay in addition to embryo morphology in embryo selection for transfer in IVF treatment if there are other embryos with the same morphology.

Furthermore, monitoring of sHLA-G in amniotic fluid and plasma of pregnant women may have an important prognostic value to recognize pathological situations.

Other interesting observations suggest that HLA-G molecules seem to be directly involved in transplant acceptance, and their analysis should be taken into consideration when monitoring transplant-patients status. In addition, soluble HLA-G plasma levels are increased in lymphoproliferative disorders or in patients suffering from malignant melanoma, glioma, breast and ovarian cancer.

Areas of investigation:

Immune Response, Infection and Inflammation

Reproduction

Transplantation

4. TEST PRINCIPLE

In the BioVendor sHLA-G ELISA, calibrators and samples are incubated in microplate wells pre-coated with monoclonal anti-sHLA-G antibody. After 60 minutes incubation and washing, monoclonal anti-human β 2-microglobulin antibody labelled with horseradish peroxidase (HRP) is added to the wells and incubated for 60 minutes with captured sHLA-G. Following another washing step, the remaining HRP conjugate is allowed to react with the substrate solution (TMB). The reaction is stopped by addition of acidic solution and absorbance of the resulting yellow product is measured. The absorbance is proportional to the concentration of sHLA-G. A calibration curve is constructed by plotting absorbance values against concentrations of calibrators, and concentrations of unknown samples are determined using this calibration curve.

5. PRECAUTIONS

- **For professional use only**
- Wear gloves and laboratory coats when handling immunodiagnostic materials
- Do not drink, eat or smoke in the areas where immunodiagnostic materials are being handled
- This kit contains components of animal origin. These materials should be handled as potentially infectious
- Avoid contact with the acidic Stop Solution and Substrate Solution, which contains hydrogen peroxide and tetramethylbenzidine (TMB). Wear gloves and eye and clothing protection when handling these reagents. Stop and/or Substrate Solutions may cause skin/eyes irritation. In case of contact with the Stop Solution and the Substrate Solution wash skin/eyes thoroughly with water and seek medical attention, when necessary
- The materials must not be pipetted by mouth

6. TECHNICAL HINTS

- Reagents with different lot numbers should not be mixed
- Use thoroughly clean glassware
- Use deionized (distilled) water, stored in clean containers
- Avoid any contamination among samples and reagents. For this purpose, disposable tips should be used for each sample and reagent
- Substrate Solution should remain colourless until added to the plate. Keep Substrate Solution protected from light
- Stop Solution should remain colourless until added to the plate. The colour developed in the wells will turn from blue to yellow immediately after the addition of the Stop Solution. Wells that are green in colour indicate that the Stop Solution has not mixed thoroughly with the Substrate Solution
- Dispose of consumable materials and unused contents in accordance with applicable national regulatory requirements

7. REAGENT SUPPLIED

<i>Kit Components</i>	<i>State</i>	<i>Quantity</i>
Antibody Coated Microtiter Strips	ready to use	96 wells
Conjugate Solution	ready to use	13 ml
Master Calibrator	lyophilized	2 vials
Dilution Buffer	ready to use	20 ml
Wash Solution Conc. (10x)	concentrated	100 ml
Substrate Solution	ready to use	13 ml
Stop Solution	ready to use	13 ml
Product Data Sheet + Certificate of Analysis	-	1 pc

8. MATERIAL REQUIRED BUT NOT SUPPLIED

- Deionized (distilled) water
- Test tubes for diluting samples
- Glassware (graduated cylinder and bottle) for Wash Solution
- Precision pipettes to deliver 10-1000 μ l with disposable tips
- Multichannel pipette to deliver 100 μ l with disposable tips
- Absorbent material (e.g. paper towels) for blotting the microtiter plate after washing
- Vortex mixer
- Orbital microplate shaker capable of approximately 300 rpm
- Microplate washer (optional). [Manual washing is possible but not preferable.]
- Microplate reader with 450 ± 10 nm filter, preferably with reference wavelength 630 nm (alternatively another one from the interval 550-650nm)
- Software package facilitating data generation and analysis (optional)

9. PREPARATION OF REAGENTS

- All reagents need to be brought to room temperature prior to use
- Always prepare only the appropriate quantity of reagents for your test
- Do not use components after the expiration date marked on their label

- Assay reagents supplied ready to use:

Antibody Coated Microtiter Strips

Stability and storage:

Return the unused strips to the provided aluminium zip-sealed bag with desiccant and seal carefully. Remaining Microtiter Strips are stable 3 months when stored at 2-8°C and protected from the moisture.

Conjugate Solution

Dilution Buffer

Substrate Solution

Stop Solution

Stability and storage:

Opened reagents are stable 3 months when stored at 2-8°C.

- Assay reagents supplied lyophilized or concentrated:

sHLA-G Master Calibrator

Refer to the Certificate of Analysis for current volume of distilled water needed for reconstitution of calibrator!!!

Reconstitute the lyophilized Master Calibrator with distilled water just prior to the assay. Let it dissolve at least 15 minutes with occasional gentle shaking (not to foam). The resulting concentration of the sHLA-G in the stock solution is **625 Units/ml**.

Prepare set of calibrators using Dilution Buffer as follows (example for duplicates):

<i>Volume of Calibrator</i>	<i>Dilution Buffer</i>	<i>Concentration</i>
Stock	-	625.00 Units/ml
100 µl of stock	400 µl	125.00 Units/ml
250 µl of 125.00 Units/ml	250 µl	62.50 Units/ml
250 µl of 62.50 Units/ml	250 µl	31.25 Units/ml
250 µl of 31.25 Units/ml	250 µl	15.63 Units/ml
250 µl of 15.63 Units/ml	250 µl	7.81 Units/ml
250 µl of 7.81 Units/ml	250 µl	3.91 Units/ml

Prepared Calibrators are ready to use, do not dilute them.

Stability and storage:

Do not store the Calibrator stock solution and set of calibrators.

Wash Solution Conc. (10x)

Dilute Wash Solution Concentrate (10x) ten-fold in distilled water to prepare a 1x working solution. Example: 100 ml of Wash Solution Concentrate (10x) + 900 ml of distilled water for use of all 96-wells.

Stability and storage:

The diluted Wash Solution is stable 1 month when stored at 2-8°C. Opened Wash Solution Concentrate (10x) is stable 3 months when stored at 2-8°C.

10. PREPARATION OF SAMPLES

The kit measures sHLA-G in serum, plasma, amniotic fluid or culture supernatant. Samples should be assayed immediately after collection or should be stored at -20°C. Mix thoroughly thawed samples just prior to the assay and avoid repeated freeze/thaw cycles, which may cause erroneous results. Avoid using hemolyzed or lipemic samples.

Serum and plasma samples:

Use undiluted serum or plasma samples.

Amniotic fluid samples:

Dilute samples 4x with Dilution Buffer just prior to the assay, e.g. 30 µl of sample + 90 µl of Dilution Buffer for singlets, or preferably 60 µl of sample + 180 µl of Dilution Buffer for duplicates. **Mix well** (not to foam). Vortex is recommended.

Stability and storage:

Samples should be stored at -20°C, or preferably at -70°C for long-term storage.

Do not store the diluted samples.

Note: It is recommended to use a precision pipette and a careful technique to perform the dilution in order to get precise results.

11. ASSAY PROCEDURE

1. Pipet **100 µl** of Calibrators, samples and Dilution Buffer (= blank), preferably in duplicates, into the appropriate wells. See *Figure 1* for example of work sheet.
2. **Incubate the plate at 2-8°C for 16-20 hours, without shaking.**
3. Wash the wells 5-times with Wash Solution (0.35 ml per well). After final wash, invert and tap the plate strongly against paper towel.
4. Add **100 µl** of Conjugate Solution into each well.
5. Incubate the plate at room temperature (ca. 25°C) for **1 hour**, shaking at ca. 300 rpm on an orbital microplate shaker.
6. Wash the wells 5-times with Wash Solution (0.35 ml per well). After final wash, invert and tap the plate strongly against paper towel.
7. Add **100 µl** of Substrate Solution into each well. Avoid exposing the microtiter plate to direct sunlight. Covering the plate with e.g. aluminium foil is recommended.
8. Incubate the plate for **25 minutes** at room temperature. The incubation time may be extended [up to 30 minutes] if the reaction temperature is below than 20°C. No shaking!
9. Stop the colour development by adding **100 µl** of Stop Solution.
10. Determine the absorbance of each well using a microplate reader set to 450 nm, preferably with the reference wavelength set to 630 nm (acceptable range: 550 – 650 nm). Subtract readings at 630 nm (550 - 650 nm) from the readings at 450 nm.
The absorbance should be read within 5 minutes following step 9.

Note: If some samples and calibrator/s have absorbances above the upper limit of your microplate reader, perform a second reading at 405 nm. A new calibrator curve, constructed using the values measured at 405 nm, is used to determine sHLA-G concentration of off-scale calibrators and samples. The readings at 405 nm should not replace the readings for samples that were “in range” at 450 nm.

Note 2: Manual washing: Aspirate wells and pipet 0.35 ml Wash Solution into each well. Aspirate wells and repeat twice. After final wash, invert and tap the plate strongly against paper towel. Make certain that Wash Solution has been removed entirely.

	strip 1+2	strip 3+4	strip 5+6	strip 7+8	strip 9+10	strip 11+12
A	Calibrator 125.00	Sample 2	Sample 10	Sample 18	Sample 26	Sample 34
B	Calibrator 62.50	Sample 3	Sample 11	Sample 19	Sample 27	Sample 35
C	Calibrator 31.25	Sample 4	Sample 12	Sample 20	Sample 28	Sample 36
D	Calibrator 15.63	Sample 5	Sample 13	Sample 21	Sample 29	Sample 37
E	Calibrator 7.81	Sample 6	Sample 14	Sample 22	Sample 30	Sample 38
F	Calibrator 3.91	Sample 7	Sample 15	Sample 23	Sample 31	Sample 39
G	Blank	Sample 8	Sample 16	Sample 24	Sample 32	Sample 40
H	Sample 1	Sample 9	Sample 17	Sample 25	Sample 33	Sample 41

Figure 1: Example of a work sheet.

12. CALCULATIONS

Most microplate readers perform automatic calculations of analyte concentration. The calibration curve is constructed by plotting the mean absorbance (Y) of Calibrators against the known concentration (X) of Calibrators in logarithmic scale, using the four-parameter algorithm. Results are reported as concentration of sHLA-G Units/ml in samples.

Alternatively, the logit log function can be used to linearize the calibration curve, i.e. logit of the mean absorbance (Y) is plotted against log of the known concentration (X) of Calibrators.

Serum or plasma samples and Calibrators are all undiluted, so there is no need to take into account any dilution factor for calculations.

The measured concentration of amniotic fluid samples calculated from the calibration curve must be multiplied by their respective dilution factor, because samples have been diluted prior to the assay, e.g. 25 U/ml (from calibration curve) x 4 (dilution factor) = 100 U/ml.

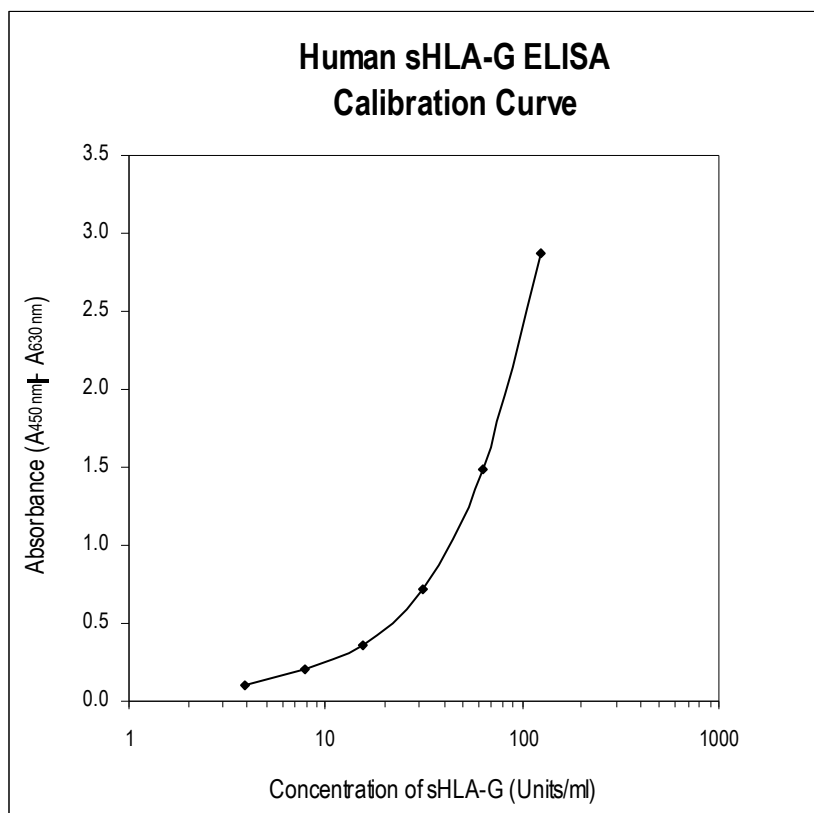


Figure 2: Typical Calibration Curve for sHLA-G ELISA.

13. PERFORMANCE CHARACTERISTICS

»» Typical analytical data of BioVendor Human sHLA-G ELISA are presented in this chapter

- **Sensitivity**

Limit of Detection (LOD), defined as concentration of analyte giving absorbance higher than mean absorbance of blank* plus three calibrator deviations of the absorbance of blank: $A_{\text{blank}} + 3 \times \text{SD}_{\text{blank}}$, is calculated from the real sHLA-G values in wells and is 3 Units/ml.

*Dilution Buffer is pipetted into blank wells.

- **Limit of assay**

Samples with absorbance exceeding the highest calibrator should be repeated with more diluted samples. Dilution factor needs to be taken into consideration in calculating the sHLA-G concentration.

- **Specificity**

Sera of several mammalian species were measured in the assay. See results below.

For details please contact us at info@biovendor.com.

<i>Mammalian serum sample</i>	<i>Observed crossreactivity</i>
Bovine	yes
Cat	no
Dog	yes
Goat	yes
Hamster	no
Horse	no
Monkey	yes
Mouse	yes
Pig	no
Rabbit	yes
Rat	no
Sheep	yes

- **Precision**

Intra-assay (Within-Run) (n=8)

<i>Sample</i>	<i>Mean (Unit/ml)</i>	<i>SD (Unit/ml)</i>	<i>CV (%)</i>
1	49.41	4.94	10.0
2	43.08	2.26	5.2

Inter-assay (Run-to-Run) (n=11)

<i>Sample</i>	<i>Mean (Unit/ml)</i>	<i>SD (Unit/ml)</i>	<i>CV (%)</i>
1	53.73	3.33	6.2
2	44.13	2.45	5.5

- **Stability of samples stored at 2-8°C**

Samples should be stored at –20°C. However, no decline in concentration of sHLA-G was observed in serum and plasma samples after 7 days when stored at 2-8°C. To avoid microbial contamination, samples were treated with ε-aminocaproic acid and sodium azide, resulting in the final concentration of 0.03% and 0.1%, respectively.

- **Effect of Freezing/Thawing**

No decline was observed in concentration of sHLA-G in serum and plasma samples after repeated (5x) freeze/thaw cycles. However it is recommended to avoid unnecessary repeated freezing/thawing of the samples.

<i>Sample</i>	<i>Number of f/t cycles</i>	<i>Serum (Unit/ml)</i>	<i>Plasma (Unit/ml)</i>		
			<i>EDTA</i>	<i>Citrate</i>	<i>Heparin</i>
1	1x	13.0	16.2	12.7	15.4
	3x	12.7	17.3	11.8	14.6
	5x	12.3	15.5	10.9	14.3
2	1x	42.2	52.2	47.1	52.1
	3x	44.6	54.6	43.8	50.5
	5x	41.3	51.3	42.5	57.8
3	1x	30.3	38.3	27.3	32.8
	3x	32.9	36.9	27.6	36.8
	5x	34.5	35.5	27.6	33.3

14. DEFINITION OF THE CALIBRATOR

The change in procedure starting the lot RD-1789R1 brought out a change in evaluation of sHLA-G levels in culture medium supernatants and amniotic fluids. Thus, the previous results (the lots up to RD-1789) should be divided by 2.0 to be comparable with the new lots starting the lot RD-1789R1.

Serum and plasma samples keeps sHLA-G value among all lots.

15. METHOD COMPARISON

The RD194070100R Human sHLA-G ELISA kit has not been compared to any other immunoassay.

16. TROUBLESHOOTING AND FAQs

»» Weak signal in all wells

Possible explanations:

- Omission of a reagent or a step
- Improper preparation or storage of a reagent
- Assay performed before reagents were allowed to come to room temperature
- Improper wavelength when reading absorbance

»» High signal and background in all wells

Possible explanations:

- Improper or inadequate washing
- Overdeveloping; incubation time with Substrate Solution should be decreased before addition of Stop Solution
- Incubation temperature over 30°C

»» High coefficient of variation (CV)

Possible explanation:

- Improper or inadequate washing
- Improper mixing Calibrators, Quality Controls or samples

17. REFERENCES

»» References to sHLA-G:

- Rebman V., Pfeiffer K., Passler M., Ferrone S., Maier S., Weiss E., Grosse-Wilde H.: Detection of soluble HLA-G molecules in plasma and amniotic fluid. *Tissue Antigens*. 1999; 53: 14-22
- Fuzzi B., Rizzo R., Criscuoli L., Noci I., Melchiorri L., Scarselli B., Bencini E., Menicucci A., Baricordi O.R.: HLA-G expression in early embryos is a fundamental prerequisite for the obtainment of pregnancy. *Eur J Immunol*. 2002; 32: 311-315
- Lila N., Amrein C., Guillemain R., Chevalier P., Latremouille C., Fabiani J.N., Dausset J., Carosella E.D., Carpentier A.: Human leukocyte antigen-G expression after heart transplantation is associated with a reduced incidence of rejection. *Circulation*. 2002; 105: 1949-1954
- Steinborn A., Rebmann V., Scharf A., Sohn C., Grosse-Wilde H.: Placental abruption is associated with decreased maternal plasma levels of soluble HLA-G. *J Clin Immunol*. 2003; 23: 307-314
- Le Bouteiller P., Pizzato N., Barakonyi A., Solier C.: HLA-G, pre-eclampsia, immunity and vascular events. *J Reprod Immunol*. 2003; 59: 219-234
- Sher G., Keskinetepe L., Nouriani M., Roussev R., Batzofin J.: Expression of sHLA-G in supernatants of individually cultured 46-h embryos: a potentially valuable indicator of 'embryo competency and IVF outcome. *Reprod Biomed Online*. 2004; 9: 74-78
- Noci I., Fuzzi B., Rizzo R., Melchiorri L., Criscuoli L., Dabizzi S., Biagiotti R., Pellegrini S., Menicucci A., Baricordi O.R.: Embryonic soluble HLA-G as a marker of developmental potential in embryos. *Hum Reprod*. 2005; 20: 138-146
- Yie S.M., Balakier H., Motamedi G., Librach C.L.: Secretion of human leukocyte antigen-G by human embryos is associated with a higher in vitro fertilization pregnancy rate. *Fertil Steril*. 2005; 83: 30-36
- Rebman V., Lemaoult J., Rouas-Friess N., Carosella E.D., Grosse-Wilde H.: Report of the Wet Workshop for Quantification of Soluble HLA-G in Essen, 2004. *Hum Immunol*. 2005; 66: 853-863
- Le Rond S., Azema C., Krawice-Radanne I., Durrbach A., Guettier C., Carosella E.D., Rouas-Freiss N.: Evidence to support the role of HLA-G5 in allograft acceptance through induction of immunosuppressive/regulatory T cells. *J Immunol*. 2006; 176: 3266-3276
- Luque J., Torres M.I., Aumente M.D., Marin J., Garcia-Jurado G., Gonzalez R., Pascual D., Guerra N., Lopez-Rubio F., Alvarez-Lopez M.R., Arizon J.M., Pena J.: Soluble HLA-G in heart transplantation: their relationship to rejection episodes and immunosuppressive therapy. *Hum Immunol*. 2006; 67: 257-263
- Carosella E.D., Favier B., Rouas-Freiss N., Moreau P., Lemaoult J.: Beyond the increasing complexity of the immunomodulatory HLA-G molecule. *Blood*. 2008; 111(10):4862-70
- Apps R., Gardner L., Moffett A.: A critical look at HLA-G. *Trends Immunol*. 2008; 29(7):313-21
- Baricordi O.R., Stignani M., Melchiorri L., Rizzo R.: HLA-G and inflammatory diseases. *Inflamm Allergy Drug Targets*. 2008; 7(2):67-74









References to this product:

- Wiendl H., Feger U., Mittelbronn M., Jack C., Schreiner B., Stadelmann C., Antel J., Brueck W., Meyermann R., Baror A., Kieseier B.C., Weller M.: Expression of the immune-tolerogenic major histocompatibility molecule HLA-G in multiple sclerosis: implications for CNS immunity. *Brain*. 2005; 128(Pt 11): 2689-2704
- Desai N., Filipovits J., Singh P., Beman M., Goldfarb J.: Assessment of soluble HLA-G levels in day 3 human embryo culture media using a new ELISA kit: Is there a relationship to embryo quality, development and implantation potential? *Fertility & Sterility*. 2005; 84, Suppl 1 (S281-S282)


- Steinborn A., Varkonyi T., Scharf A., Bahlmann F., Klee A., Sohn C.: Early detection of decreased soluble HLA-G levels in the maternal circulation predicts the occurrence of preeclampsia and intrauterine growth retardation during further course of pregnancy. *Am J Reprod Immunol* . 2007; 7(4):277-86
- Tahan F., Patiroglu T. Plasma soluble human leukocyte antigen G levels in asthmatic children. *Int Arch Allergy Immunol* . 2006; 141(3):213-6 (2006)
- Desai N., Filipovits J., Goldfarb J.: Secretion of soluble HLA-G by day 3 human embryos associated with higher pregnancy and implantation rates: assay of culture media using a new ELISA kit. *Repro Biomed Online* . 2006; 13(2):272-7
- Coulam C.B., Roussev R.G.: Soluble HLA G concentrations in culture media from in vitro fertilized human embryos are predictive of implantation potential. *American Journal of Reproductive Immunology*. 2005; 53:279-309
- Wang F., Wen Z., Li H., Yang Z., Zhao X., Yao X.: Human leukocyte antigen-G is expressed by the eutopic and ectopic endometrium of adenomyosis. *Fertil Steril*. 2008; 90(5):1599-604
- Shaikly V.R., Morrison I.E., Taranissi M., Noble C.V., Withey A.D., Cherry R.J., Blois S.M. and Fernández N.: Analysis of HLA-G in maternal plasma, follicular fluid, and preimplantation embryos reveal an asymmetric pattern of expression. *J Immunol*. 2008; 180(6):4330-7
- Chen X.Y., Yan W.H., Lin A., Xu H.H., Zhang J.G., Wang X.X.: The 14 bp deletion polymorphisms in HLA-G gene play an important role in the expression of soluble HLA-G in plasma. *Tissue Antigens*. 2008; 72(4):335-41
- Ciprandi G., Colombo B.M., Contini P., Cagnati P., Pistorio A., Puppo F., Murdaca G.: Soluble HLA-G and HLA-A,-B,-C serum levels in patients with allergic rhinitis. *Allergy*. 2008; 63(10):1335-8
- Eidukaite A., Tamosiunas V.: Soluble HLA-G in the peritoneal fluid of women with endometriosis. *Fertil Steril*. 2008; 89(2):465-7
- Le Maux A., Noël G., Birebent B., Grosset J.M., Vu N., De Guibert S., Bernard M., Semana G., Amiot L.: Soluble human leukocyte antigen-G molecules in peripheral blood haematopoietic stem cell transplantation: a specific role to prevent acute graft-versus-host disease and a link with regulatory T cells. *Clin Exp Immunol*. 2008; 152(1):50-6
- Giannopoulos K., Schmitt M., Kowal M., Wlasiuk P., Bojarska-Junak A., Roliński J., Dmoszyńska A.: The significance of soluble HLA-G plasma levels as well as messenger HLA-G for B-cell chronic lymphocytic leukemia (B-CLL). *Leuk Res*. 2008; 32(12):1815-9
- McCormick J., Whitley G.S., Le Bouteiller P., Cartwright J.E.: Soluble HLA-G regulates motility and invasion of the trophoblast-derived cell line SGHPL-4. *Hum Reprod*. 2009; 24(6):1339-45
- Kusanovic J.P., Romero R., Jodicke C., Mazaki-Tovi S., Vaisbuch E., Erez O., Mittal P., Gotsch F., Chaiworapongsa T., Edwin S.S., Pacora P., Hassan S.S.: Amniotic fluid soluble human leukocyte antigen-G in term and preterm parturition, and intra-amniotic infection/inflammation. *J Matern Fetal Neonatal Med*. 2009; 22(12):1151-66
- Roussev R.G., Acacio B., Parrett S., Romineb H., Coulam C.B.: Soluble HLA-G as potential candidate to help morphology grading score in embryo selection for IVF/ET. *Journal of Reproductive Immunology*. 2009; 81, 113–175
- Ivanova-Todorova E., Mourdjeva M., Kyurkchiev D., Bochev I., Stoyanova E., Dimitrov R., Timeva T., Yunakova M., Bukarev D., Shterev A., Tivchev P., Kyurkchiev S.: HLA-G expression is up-regulated by progesterone in mesenchymal stem cells. *Am J Reprod Immunol*. 2009; 62(1):25-33
- Mach P., Blecharz P., Basta P., Marianowski P., Skret-Magierlo J., Kojs Z., Grabiec M., Wicherek L.: Differences in the soluble HLA-G blood serum concentration levels in patients with ovarian cancer and ovarian and deep endometriosis. *Am J Reprod Immunol*. 2010; 63(5):387-95
- Chen C.H., Liao H.T., Chen H.A., Liu C.H., Liang T.H., Wang C.T., Tsai C.Y., Chou C.T.: Human leukocyte antigen-G in ankylosing spondylitis and the response after tumour necrosis factor-alpha blocker therapy. *Rheumatology (Oxford)*. 2010; 49(2):264-70

»» For more references on this product see our WebPages at www.biovendor.com

18. EXPLANATION OF SYMBOLS

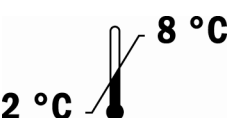	Catalogue number
	Content
	Lot number
	See instructions for use
	Biological hazard
	Expiry date
	Storage conditions
	Identification of packaging materials

Assay Procedure Summary

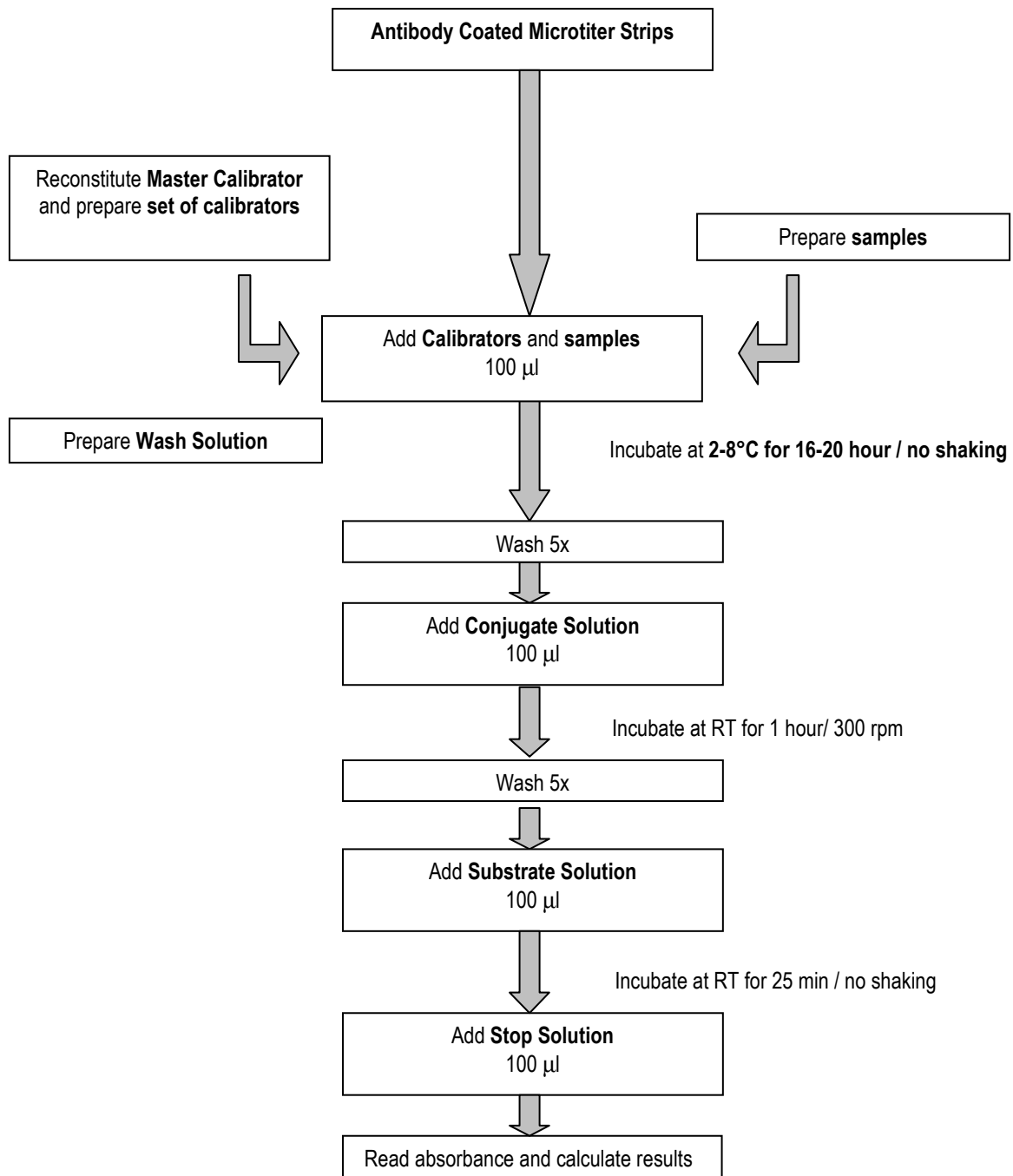

1							
2							
3							
4							
5							
6							
7							
8							
9							
10							
11							
12							
A							
B							
C							
D							
E							
F							
G							
H							

NOTES





EXBIO Praha a.s.	Nad Safinou II 366	252 42 Vestec CZECH REPUBLIC	Phone: +420-261090664 Fax: +420-261090661	E-mail: exbio@exbio.cz Web: http://www.exbio.cz
-------------------------	---------------------------	-----------------------------------------	------------------------------------------------------	----------------------------------------------------------------------------------------------------------------------------------



HEADQUARTERS: BioVendor - Laboratorní medicína a.s.	Karasek 1767/1	621 00 Brno CZECH REPUBLIC	Phone: +420-549-124-185 Fax: +420-549-211-460	E-mail: info@biovendor.com Web: www.biovendor.com
EUROPEAN UNION: BioVendor GmbH	Im Neuenheimer Feld 583	D-69120 Heidelberg GERMANY	Phone: +49-6221-433-9100 Fax: +49-6221-433-9111	E-mail: infoEU@biovendor.com
USA, CANADA AND MEXICO: BioVendor LLC	1463 Sand Hill Road Suite 227	Candler, NC 28715 USA	Phone: +1-828-670-7807 +1-800-404-7807 Fax: +1-828-670-7809	E-mail: infoUSA@biovendor.com
CHINA - Hong Kong Office: BioVendor Laboratories Ltd	Room 4008 Hong Kong Plaza, No.188	Connaught Road West Hong Kong, CHINA	Phone: +852-2803-0523 Fax: +852-2803-0525	E-mail: infoHK@biovendor.com
CHINA – Mainland Office: BioVendor Laboratories Ltd	Room 2917, 29F R & F Ying Feng Plaza, No.2 Huaqiang road	Pearl River New Town Guang Zhou, CHINA	Phone: +86-20-38065519 Fax: +86-20-38065529	E-mail: infoCN@biovendor.com