





**RUO** in the USA

Please use only the valid version of the package insert provided with the kit.

This kit is intended for Research Use Only.

Not for use in diagnostic procedures.

#### **INTENDED USE**

Competitive immunoenzymatic colorimetric method for determination of Free Cortisol concentration in Urine. Urinary "Free" Cortisol kit is intended for laboratory use only.

#### PRINCIPLE

Urinary Cortisol (antigen) in the sample competes with horseradish-peroxidase Cortisol (enzyme-labelled-antigen) for binding onto the limited number of anti-Cortisol (antibody) sites on the microplates (solid phase). After incubation, the bound/free separation is performed by a simple solid-phase washing. The enzyme substrate ( $H_2O_2$ ) and the TMB-Sustrate (TMB) are added. After an appropriate time has elapsed for maximum colour development, the enzyme reaction is stopped and the absorbances are determined. Urinary Cortisol concentration in the sample is calculated based on a series of standard. The colour intensity is inversely proportional to the urinary cortisol concentration in the sample.

#### ReagentS, materialS and instrumentation

#### Reagents and materials supplied in the kit

- 1. Cortisol Standards STD0 STD4, 5 vials, 1 mL each
- 2. **Incubation buffer,** 1 vial, 100 mL Phosphate buffer 50 mM pH 7.4; BSA 1 g/L
- 3. **Conjugate**, 1 vial, 1.0 mL Cortisol-HRP conjugate
- 4. **Coated Microplate**, 1 breakable microplate Anti-Cortisol-IgG adsorbed on microplate
- TMB-substrate, 1 vial, 15 mL H<sub>2</sub>O<sub>2</sub>-TMB 0.26 g/L (avoid any skin contact)
- 6. **Stop solution,** 1 vial, 15 mL Sulphuric acid 0.15 mol/L (avoid any skin contact)
- 7. **Low Control,** 1 vial, 1 mL Ready to use





# **C E** Revised 26 Jan. 2011 rm (Vers. 8.1)

8. **High Control,** 1 vial, 1 mL Ready to use

#### **Reagents necessary not supplied**

Distilled water

#### Auxiliary materials and instrumentation

Automatic dispenser Microplates reader (450 nm)

#### Note

Store all reagents at  $2 \degree C - 8 \degree C$  in the dark.

Open the bag of reagent 4 (Coated Microplate) only when it is at room temperature and close immediately after use. Do not remove the adhesive sheet from the unused strips.

#### PRECAUTIONS

The reagents contain Proclin 300 as preservative.

Avoid the exposure of reagent TMB/H<sub>2</sub>O<sub>2</sub> to directed sunlight, metals or oxidants.

Maximum precision is required for reconstitution and dispensation of the reagents.

Do not use different lots of reagents.

Do not use heavily haemolized samples.

This method allows the determination of Cortisol from10 ng/mL to 500 ng/mL.

The clinical significance of the Cortisol determination can be invalidated if the sample was treated with corticosteroids or natural or synthetic steroids.



in the USA





## **C E** Revised 26 Jan. 2011 rm (Vers. 8.1)

**RUO** in the USA

### PROCEDURE

#### Preparation of the Standard (S<sub>0</sub>, S<sub>1</sub>, S<sub>2</sub>, S<sub>3</sub>, S<sub>4</sub>)

The standards have the following concentration of Cortisol:							
	$\mathbf{S}_{0}$	$\mathbf{S}_1$	$S_2$	S <sub>3</sub>	$S_4$		
ng/mL	0	10	50	150	500		

	0				
C	table when	stard at $\perp 1$	°C until the	avniration	data of

Stable when stored at +4 °C until the expiration date of the kit; once opened the standards are stable six months at 2 °C - 8 °C.

#### Preparation of diluted Conjugate

Prepare immediately before use.

Add 10 µL Conjugate (reagent 3) to 1.0 mL of Incubation Buffer (reagent 2). Mix gently for at least 10 minutes with a rotating mixer.

Stable for 3 hours at room temperature (22 °C – 28 °C).

#### Preparation of the Sample

The total volume of urine excreted during a 24 hours should be collected and mixed in a single container. Urine samples which are not to be assayed immediately should be stored at  $2 \degree C - 8 \degree C$  or at  $-20\degree C$  for periods longer than a week.

Dilute the urine sample (1 + 1) with Incubation buffer (e.g. 100  $\mu$ L + 100  $\mu$ L)





## **C E** Revised 26 Jan. 2011 rm (Vers. 8.1)

## **RUO** in the USA

#### Procedure

As it is necessary to perform the determination in duplicate, prepare two wells for each of the four points of the standard curve  $(S_0-S_4)$ , one for Blank.

Reagent	Standards/Controls	Samples	Blank			
Standard S0-S4 Controls	10 µL					
Diluted samples		10 μL				
Diluted Conjugate	300 μL	300 μL				
Incubate at 37°C for 1 hour. Remove the contents from each well; wash the wells 2 times with 400 μL of distilled water.						
TMB substrate	100 μL	100 µL	100 µL			
Incubate at room temperature (22 °C – 28 °C) for 15 minutes in the dark						
Stop solution	100 μL	100 µL	100 µL			
Shake the microplate gently. Read the absorbance (E) at 450 nm against Blank.						

### QUALITY CONTROL

Each laboratory should assay controls at normal, high and low levels range of Urinary Cortisol for monitoring assay performance. These controls should be treated as unknowns and values determined in every test procedure performed. Quality control charts should be maintained to follow the performance of the supplied reagents. Pertinent statistical methods should be employed to ascertain trends. The individual laboratory should set acceptable assay performance limits. Other parameters that should be monitored include the 80, 50 and 20% intercepts of the standard curve for run-to-run reproducibility. In addition, maximum absorbance should be consistent with past experience. Significant deviation from established performance can indicate unnoticed change in experimental conditions or degradation of kit reagents. Fresh reagents should be used to determine the reason for the variations.

### LIMITATIONS OF PROCEDURE

#### **Assay Performance**

Sample(s), which are contaminated microbiologically, should not be used in the assay.

Highly lipemeic or haemolysed specimen(s) should similarly not be used.

It is important that the time of reaction in each well is held constant for reproducible results. Pipetting of samples should not extend beyond ten minutes to avoid assay drift. If more than one plate is used, it is recommended to repeat the dose response curve.





Addition of the substrate solution initiates a kinetic reaction, which is terminated by the addition of the stop solution. Therefore, the addition of the substrate and the stopping solution should be added in the same sequence to eliminate any time deviation during reaction.

Plate readers measure vertically. Do not touch the bottom of the wells.

Failure to remove adhering solution adequately in the aspiration or decantation wash step(s) may result in poor replication and spurious results.

#### Interpretation

If computer controlled data reduction is used to calculate the results of the test, it is imperative that the predicted values for the calibrators fall within 10% of the assigned concentrations.

#### RESULTS

#### **Mean Absorbance**

Calculate the mean of the absorbance (Em) for each point of the standard curve and of each sample

#### **Standard Curve**

Plot the values of absorbance of the standards against concentration. Draw the best-fit curve through the plotted points (e.g.: Four Parameter Logistic).

#### **Calculation of Results**

Interpolate the values of the samples on the standard curve to obtain the corresponding values of the concentrations expressed in ng/mL.

To calculate the cortisol concentration in urine, calculate as above and correct for total volume of volume of urine collected in 24 hours:

ng/mL x Vol(mL) urine 24 h/  $1.000 = \mu g$  Cortisol/24h

#### **Reference Values**

 $50 - 190 \ \mu g/24 \ hours$ 



**RUO** in the USA





# **C E** Revised 26 Jan. 2011 rm (Vers. 8.1)

**RUO** in the USA

#### WASTE MANAGEMENT

Reagents must be disposed off in accordance with local regulations.

#### BIBLIOGRAPHY

- 1. Foster, L. B. and Dunn, R. T. Clin. Chem. 20/3 365(1974)
- 2. De Lacerda, et al J. Clin. Endocr. and Metab. 36,227 (1973)
- 3. Rolleri, E., et al Clin. Chim. Acta 66 319 (1976)
- 4. Kobayashi, Y., et al Steroids, 32 no 1(1978)
- 5. Akarawa, et al Anal. Biochem. 97 248 (1979)





# **C E** Revised 26 Jan. 2011 rm (Vers. 8.1)

RUO in

in the USA

#### TROUBLESHOOTING

#### ERRORS / POSSIBLE CAUSES / SUGGESTIONS

- No colorimetric reaction
- no conjugate pipetted
- contamination of conjugates and/or of substrate
- errors in performing the assay procedure (e.g. accidental pipetting of reagents in a wrong sequence or from the wrong vial, etc.)

#### Too low reaction (too low ODs)

- incorrect conjugate (e.g. not from original kit)
- incubation time too short, incubation temperature too low

#### Too high reaction (too high ODs)

- incorrect conjugate (e.g. not from original kit)
- incubation time too long, incubation temperature too high
- water quality for wash buffer insufficient (low grade of deionization)
- insufficient washing (conjugates not properly removed)

#### **Unexplainable outliers**

- contamination of pipettes, tips or containers
- insufficient washing (conjugates not properly removed)
- too high within-run CV%
- reagents and/or strips not pre-warmed to room temperature prior to use
- plate washer is not washing correctly (suggestion: clean washer head)
- too high between-run CV %
- incubation conditions not constant (time, temperature)
- controls and samples not dispensed at the same time (with the same intervals) (check pipetting order)
- person-related variation

#### BIBLIOGRAPHY

- 1. Foster, L.B. and Dunn, R.T. Clin Chem: 20/3, 365 (1974)
- 2. De Laceda, L., Kowarski, A., and Migeon, C.J. J. Clin Endocr. and Metab: 36:227 (1973)
- 3. Rolleri, E., Zannino, M., Orlandini, S., Malvano, R. Clin chim Acta 66 319 (1976)
- 4. Kobayashi, Y., et al. Steroids, 32 no. 1 (1978)
- 5. Arakawa, H., Maeda, M., Tsuji, A. Anal. Biochem. 97248 (1979)