

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

Please use only the valid version of the package insert provided with the kit.

This kit is intended for Research Use Only.

This kit is not intended for in vitro diagnostic procedures.

INTENDED USE

The pig IFN- γ ELISA is an enzyme-linked immunosorbent assay for detection of pig IFN- γ .

The pig IFN- γ ELISA is for research use only. Not for diagnostic or therapeutic procedures.

PRINCIPLES OF THE TEST

An anti-pig IFN- γ coating antibody is adsorbed onto microwells.

Pig IFN- γ present in the sample or standard binds to antibodies adsorbed to the microwells.

Following incubation unbound biological components are removed during a wash step. A biotin-conjugated anti-pig IFN- γ antibody is added and binds to pig IFN- γ captured by the first antibody.

Following incubation unbound biotin-conjugated anti-pig IFN- γ antibody is removed during a wash step. Streptavidin-HRP is added and binds to the biotin-conjugated anti-pig IFN- γ antibody.

Following incubation unbound Streptavidin-HRP is removed during a wash step, and substrate solution reactive with HRP is added to the wells.

A coloured product is formed in proportion to the amount of pig IFN- γ present in the sample or standard. The reaction is terminated by addition of acid and absorbance is measured at 450 nm. A standard curve is prepared from 7 pig IFN- γ standard dilutions and pig IFN- γ sample concentration determined.

REAGENTS PROVIDED

Reagents for pig IFN- γ ELISA (96 tests)

- 1 aluminium pouch with a **Microwell Plate coated** with polyclonal antibody to pig IFN- γ
- 1 vial (150 μ l) **Biotin-Conjugate** anti-pig IFN- γ monoclonal antibody
- 1 vial (150 μ l) **Streptavidin-HRP**
- 2 vials pig IFN- γ **Standard** lyophilized, 4000 pg/ml upon reconstitution
- 1 vial (5 ml) **Assay Buffer Concentrate** 20x (PBS with 1% Tween 20 and 10% BSA)
- 1 bottle (50 ml) **Wash Buffer Concentrate** 20x (PBS with 1% Tween 20)
- 1 vial (15 ml) **Substrate Solution** (tetramethyl-benzidine)
- 1 vial (12 ml) **Stop Solution** (1M Phosphoric acid)

DRG[®] Interferon γ (porcine) (EIA-4828)**Revised 27 Jan. 2011 rm (Vers. 3.1)****For Veterinary Use Only**

- 1 vial (0.4 ml) **Blue-Dye**
- 1 vial (0.4 ml) **Green-Dye**
- 1 vial (0.4 ml) **Red-Dye**
- 6 **Adhesive Films**

STORAGE INSTRUCTIONS – ELISA KIT

Store kit reagents between 2°C and 8°C. Immediately after use remaining reagents should be returned to cold storage (2°C to 8°C). Expiry of the kit and reagents is stated on labels.

Expiry of the kit components can only be guaranteed if the components are stored properly, and if, in case of repeated use of one component, this reagent is not contaminated by the first handling.

SPECIMEN COLLECTION AND STORAGE INSTRUCTIONS

Cell culture supernatant and serum were tested with this assay. Other biological samples might be suitable for use in the assay. Remove serum from the clot as soon as possible after clotting.

Samples containing a visible precipitate must be clarified prior to use in the assay. Do not use grossly hemolyzed or lipemic specimens.

Samples should be aliquoted and must be stored frozen at -20°C to avoid loss of bioactive pig IFN- γ . If samples are to be run within 24 hours, they may be stored at 2°C to 8°C (for sample stability refer to 0).

Avoid repeated freeze-thaw cycles. Prior to assay, the frozen sample should be brought to room temperature slowly and mixed gently.

MATERIALS REQUIRED BUT NOT PROVIDED

- 5 ml and 10 ml graduated pipettes
- 5 μ l to 1000 μ l adjustable single channel micropipettes with disposable tips
- 50 μ l to 300 μ l adjustable multichannel micropipette with disposable tips
- Multichannel micropipette reservoir
- Beakers, flasks, cylinders necessary for preparation of reagents
- Device for delivery of wash solution (multichannel wash bottle or automatic wash system)
- Microwell strip reader capable of reading at 450 nm (620 nm as optional reference wave length)
- Glass-distilled or deionized water
- Statistical calculator with program to perform regression analysis

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only**PRECAUTIONS FOR USE**

- All chemicals should be considered as potentially hazardous. We therefore recommend that this product is handled only by those persons who have been trained in laboratory techniques and that it is used in accordance with the principles of good laboratory practice. Wear suitable protective clothing such as laboratory overalls, safety glasses and gloves. Care should be taken to avoid contact with skin or eyes. In the case of contact with skin or eyes wash immediately with water. See material safety data sheet(s) and/or safety statement(s) for specific advice.
- Reagents are intended for research use only and are not for use in diagnostic or therapeutic procedures.
- Do not mix or substitute reagents with those from other lots or other sources.
- Do not use kit reagents beyond expiration date on label.
- Do not expose kit reagents to strong light during storage or incubation.
- Do not pipette by mouth.
- Do not eat or smoke in areas where kit reagents or samples are handled.
- Avoid contact of skin or mucous membranes with kit reagents or specimens.
- Rubber or disposable latex gloves should be worn while handling kit reagents or specimens.
- Avoid contact of substrate solution with oxidizing agents and metal.
- Avoid splashing or generation of aerosols.
- In order to avoid microbial contamination or cross-contamination of reagents or specimens which may invalidate the test use disposable pipette tips and/or pipettes.
- Use clean, dedicated reagent trays for dispensing the conjugate and substrate reagent.
- Exposure to acid inactivates the conjugate.
- Glass-distilled water or deionized water must be used for reagent preparation.
- Substrate solution must be at room temperature prior to use.
- Decontaminate and dispose specimens and all potentially contaminated materials as they could contain infectious agents. The preferred method of decontamination is autoclaving for a minimum of 1 hour at 121.5°C.
- Liquid wastes not containing acid and neutralized waste may be mixed with sodium hypochlorite in volumes such that the final mixture contains 1.0% sodium hypochlorite. Allow 30 minutes for effective decontamination. Liquid waste containing acid must be neutralized prior to the addition of sodium hypochlorite.

PREPARATION OF REAGENTS

Buffer Concentrates should be brought to room temperature and should be diluted before starting the test procedure. If crystals have formed in the **Buffer Concentrates**, warm them gently until they have completely dissolved.

Wash Buffer (1x)

Pour entire contents (50 ml) of the **Wash Buffer Concentrate** (20x) into a clean 1000 ml graduated cylinder. Bring to final volume of 1000 ml with glass-distilled or deionized water. Mix gently to avoid foaming. The pH of the final solution should adjust to 7.4.

Transfer to a clean wash bottle and store at 2°C to 25°C. Please note that Wash Buffer (1x) is stable for 30 days.

Wash Buffer (1x) may also be prepared as needed according to the following table:

DRG International Inc., USA

Fax: (908) 233-0758 • **E-mail:** corp@drg-international.com • **Web:** www.drg-international.com

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

Number of Strips	Wash Buffer Concentrate (20x) (ml)	Distilled Water (ml)
1 - 6	25	475
1 - 12	50	950

Assay Buffer (1x)

Pour the entire contents (5 ml) of the **Assay Buffer Concentrate** (20x) into a clean 100 ml graduated cylinder. Bring to final volume of 100 ml with distilled water. Mix gently to avoid foaming.

Store at 2°C to 8°C. Please note that the Assay Buffer (1x) is stable for 30 days.

Assay Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Assay Buffer Concentrate (20x) (ml)	Distilled Water (ml)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

Biotin-Conjugate

Please note that the Biotin-Conjugate should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated **Biotin-Conjugate** solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Biotin-Conjugate (ml)	Assay Buffer (1x) (ml)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

Streptavidin-HRP

Please note that the Streptavidin-HRP should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated **Streptavidin-HRP** solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Streptavidin-HRP (ml)	Assay Buffer (1x) (ml)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

Pig IFN- γ Standard

Reconstitute **pig IFN- γ standard** by addition of distilled water.

Reconstitution volume is stated on the label of the standard vial. Swirl or mix gently to insure complete and homogeneous solubilization (concentration of reconstituted standard = 4000 pg/ml).

Allow the reconstituted standard to sit for a minimum of 10 minutes. Mix well prior to making dilutions.

After usage remaining standard cannot be stored and has to be discarded.

Standard dilutions can be prepared directly on the microwell plate (see 10.c) or alternatively in tubes (see 1.1.1).

1.1.1 External Standard Dilution

Label 7 tubes, one for each standard point.

S1, S2, S3, S4, S5, S6, S7

Then prepare 1:2 serial dilutions for the standard curve as follows:

Pipette 225 μ l of Assay Buffer (1x) into each tube.


Pipette 225 μ l of reconstituted standard (concentration = 4000 pg/ml) into the first tube, labelled S1, and mix (concentration of standard 1 = 2000 pg/ml).

Pipette 225 μ l of this dilution into the second tube, labelled S2, and mix thoroughly before the next transfer.

Repeat serial dilutions 5 more times thus creating the points of the standard curve (see Figure 1).

Assay Buffer (1x) serves as blank.

Figure 1

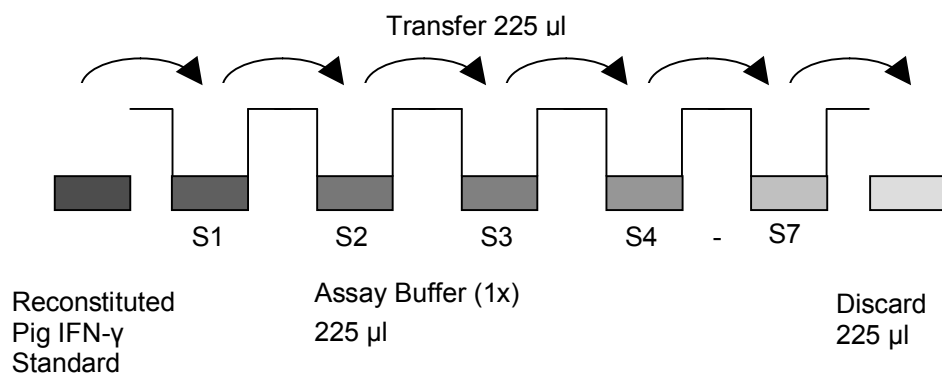

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only**Addition of Colour-giving Reagents: Blue-Dye, Green-Dye, Red-Dye**

In order to help our customers to avoid any mistakes in pipetting the ELISAs, a tool is offered that helps to monitor the addition of even very small volumes of a solution to the reaction well by giving distinctive colours to each step of the ELISA procedure.

This procedure is optional, does not in any way interfere with the test results, and is designed to help the customer with the performance of the test, but can also be omitted, just following the instruction booklet.

Alternatively, the dye solutions from the stocks provided (*Blue-Dye*, *Green-Dye*, *Red-Dye*) can be added to the reagents according to the following guidelines:

- 1. Diluent:** Before standard and sample dilution add the *Blue-Dye* at a dilution of 1:250 (see table below) to the appropriate diluent (1x) according to the test protocol. After addition of *Blue-Dye*, proceed according to the instruction booklet.

5 ml Assay Buffer (1x)	20 μ l <i>Blue-Dye</i>
12 ml Assay Buffer (1x)	48 μ l <i>Blue-Dye</i>
50 ml Assay Buffer (1x)	200 μ l <i>Blue-Dye</i>

- 2. Biotin-Conjugate:** Before dilution of the concentrated Biotin-Conjugate, add the *Green-Dye* at a dilution of 1:100 (see table below) to the Assay Buffer (1x) used for the final conjugate dilution. Proceed after addition of *Green-Dye* according to the instruction booklet: Preparation of Biotin-Conjugate.

6 ml Assay Buffer (1x)	60 μ l <i>Green-Dye</i>
12 ml Assay Buffer (1x)	120 μ l <i>Green-Dye</i>

- 3. Streptavidin-HRP:** Before dilution of the concentrated Streptavidin-HRP, add the *Red-Dye* at a dilution of 1:250 (see table below) to the Assay Buffer (1x) used for the final Streptavidin-HRP dilution. Proceed after addition of *Red-Dye* according to the instruction booklet: Preparation of Streptavidin-HRP.

6 ml Assay Buffer (1x)	24 μ l <i>Red-Dye</i>
12 ml Assay Buffer (1x)	48 μ l <i>Red-Dye</i>

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

TEST PROTOCOL

- Determine the number of microwell strips required to test the desired number of samples plus appropriate number of wells needed for running blanks and standards. Each sample, standard, blank and optional control sample should be assayed in duplicate. Remove extra microwell strips from holder and store in foil bag with the desiccant provided at 2°C-8°C sealed tightly.
- Wash the microwell strips twice with approximately 400 μ l **Wash Buffer** per well with thorough aspiration of microwell contents between washes. Allow the Wash Buffer to sit in the wells for about **10 – 15 seconds** before aspiration. Take care not to scratch the surface of the microwells.
After the last wash step, empty wells and tap microwell strips on absorbent pad or paper towel to remove excess Wash Buffer. Use the microwell strips immediately after washing. Alternatively microwell strips can be placed upside down on a wet absorbent paper for not longer than 15 minutes. **Do not allow wells to dry.**
- Standard dilution on the microwell plate** (Alternatively the standard dilution can be prepared in tubes - see 1.1.1):
Add 100 μ l of Assay Buffer (1x) in duplicate to all **standard wells**. Pipette 100 μ l of prepared **standard** (see Preparation of Standard 0, concentration = 4000 pg/ml) in duplicate into well A1 and A2 (see Table 1). Mix the contents of wells A1 and A2 by repeated aspiration and ejection (concentration of standard 1, S1 = 2000 pg/ml), and transfer 100 μ l to wells B1 and B2, respectively (see Figure 2). Take care not to scratch the inner surface of the microwells. Continue this procedure 5 times, creating two rows of pig IFN- γ standard dilutions ranging from 2000 to 31 pg/ml. Discard 100 μ l of the contents from the last microwells (G1, G2) used.

Figure 2

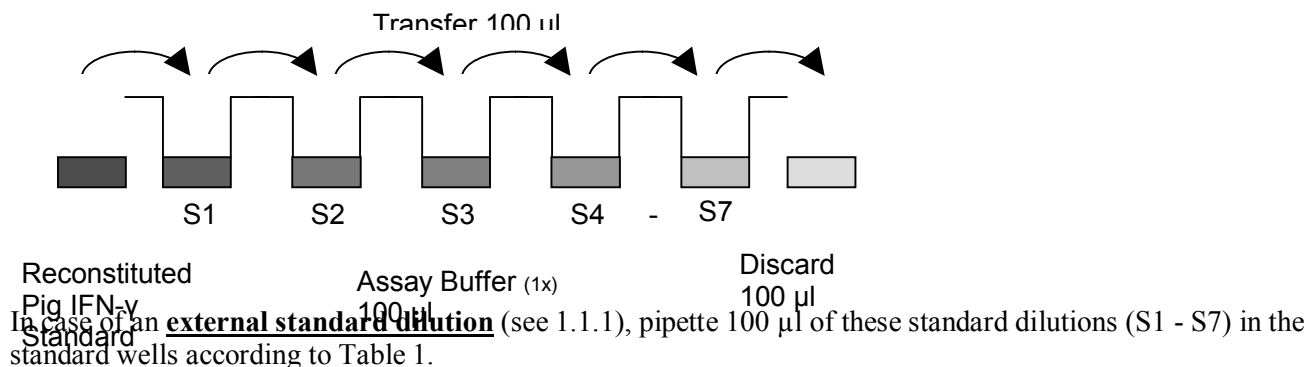


Table 1

Table depicting an example of the arrangement of blanks, standards and samples in the microwell strips:

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

	1	2	3	4
A	Standard 1 (2000 pg/ml)	Standard 1 (2000 pg/ml)	Sample 1	Sample 1
B	Standard 2 (1000 pg/ml)	Standard 2 (1000 pg/ml)	Sample 2	Sample 2
C	Standard 3 (500 pg/ml)	Standard 3 (500 pg/ml)	Sample 3	Sample 3
D	Standard 4 (250 pg/ml)	Standard 4 (250 pg/ml)	Sample 4	Sample 4
E	Standard 5 (125 pg/ml)	Standard 5 (125 pg/ml)	Sample 5	Sample 5
F	Standard 6 (63 pg/ml)	Standard 6 (63 pg/ml)	Sample 6	Sample 6
G	Standard 7 (31 pg/ml)	Standard 7 (31 pg/ml)	Sample 7	Sample 7
H	Blank	Blank	Sample 8	Sample 8

- d. Add 100 μ l of **Assay Buffer (1x)** in duplicate to the **blank wells**.
- e. Add 50 μ l of **Assay Buffer (1x)** to the **sample wells**.
- f. Add 50 μ l of each **sample** in duplicate to the **sample wells**.
- g. Cover with an adhesive film and incubate at room temperature (18 to 25°C) for 1 hour, if available on a microplate shaker set at 200 rpm.
- h. Prepare **Biotin-Conjugate** (see Preparation of Biotin-Conjugate 0).
- i. Remove adhesive film and empty wells. **Wash** microwell strips 3 times according to point b. of the test protocol. Proceed immediately to the next step.
- j. Add 100 μ l of **Biotin-Conjugate** to all wells.
- k. Cover with an adhesive film and incubate at room temperature (18 to 25°C) for 1 hour, if available on a microplate shaker set at 200 rpm.
- l. Prepare **Streptavidin-HRP** (refer to Preparation of Streptavidin-HRP 0).
- m. Remove adhesive film and empty wells. **Wash** microwell strips 3 times according to point b. of the test protocol. Proceed immediately to the next step.
- n. Add 100 μ l of diluted **Streptavidin-HRP** to all wells, including the blank wells.
- o. Cover with an adhesive film and incubate at room temperature (18° to 25°C) for 30 minutes, if available on a microplate shaker set at 200 rpm.
- p. Remove adhesive film and empty wells. **Wash** microwell strips 3 times according to point b. of the test protocol. Proceed immediately to the next step.
- q. Pipette 100 μ l of **TMB Substrate Solution** to all wells.
- r. Incubate the microwell strips at room temperature (18° to 25°C) for about 10 min. Avoid direct exposure to intense light.
The colour development on the plate should be monitored and the substrate reaction stopped (see next point of this protocol) before positive wells are no longer properly recordable.
Determination of the ideal time period for colour development has to be done individually for each assay.
 It is recommended to add the stop solution when the highest standard has developed a dark blue colour. Alternatively the colour development can be monitored by the ELISA reader at 620 nm. The substrate reaction should be stopped as soon as Standard 1 has reached an OD of 0.6 – 0.65.

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

- s. Stop the enzyme reaction by quickly pipetting 100 μ l of **Stop Solution** into each well. It is important that the Stop Solution is spread quickly and uniformly throughout the microwells to completely inactivate the enzyme. Results must be read immediately after the Stop Solution is added or within one hour if the microwell strips are stored at 2 - 8°C in the dark.
- t. Read absorbance of each microwell on a spectro-photometer using 450 nm as the primary wave length (optionally 620 nm as the reference wave length; 610 nm to 650 nm is acceptable). Blank the plate reader according to the manufacturer's instructions by using the blank wells. Determine the absorbance of both the samples and the standards.

Note: In case of incubation without shaking the obtained O.D. values may be lower than indicated below. Nevertheless the results are still valid.

CALCULATION OF RESULTS

- Calculate the average absorbance values for each set of duplicate standards and samples. Duplicates should be within 20 per cent of the mean value.
- Create a standard curve by plotting the mean absorbance for each standard concentration on the ordinate against the pig IFN- γ concentration on the abscissa. Draw a best fit curve through the points of the graph (a 5-parameter curve fit is recommended).
- To determine the concentration of circulating pig IFN- γ for each sample, first find the mean absorbance value on the ordinate and extend a horizontal line to the standard curve. At the point of intersection, extend a vertical line to the abscissa and read the corresponding pig IFN- γ concentration.
- **If instructions in this protocol have been followed samples have been diluted 1:2 (50 μ l sample + 50 μ l Assay Buffer (1x)), the concentration read from the standard curve must be multiplied by the dilution factor (x 2).**
- **Calculation of samples with a concentration exceeding standard 1 may result in incorrect pig IFN- γ levels. Such samples require further external predilution according to expected pig IFN- γ values with Assay Buffer (1x) in order to precisely quantitate the actual pig IFN- γ level.**
- It is suggested that each testing facility establishes a control sample of known pig IFN- γ concentration and runs this additional control with each assay. If the values obtained are not within the expected range of the control, the assay results may be invalid.
- A representative standard curve is shown in Figure 3. This curve cannot be used to derive test results. Each laboratory must prepare a standard curve for each group of microwell strips assayed.


Figure 3

Representative standard curve for pig IFN- γ ELISA. Pig IFN- γ was diluted in serial 2-fold steps in Assay Buffer (1x). Do not use this standard curve to derive test results. A standard curve must be run for each group of microwell strips assayed.

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

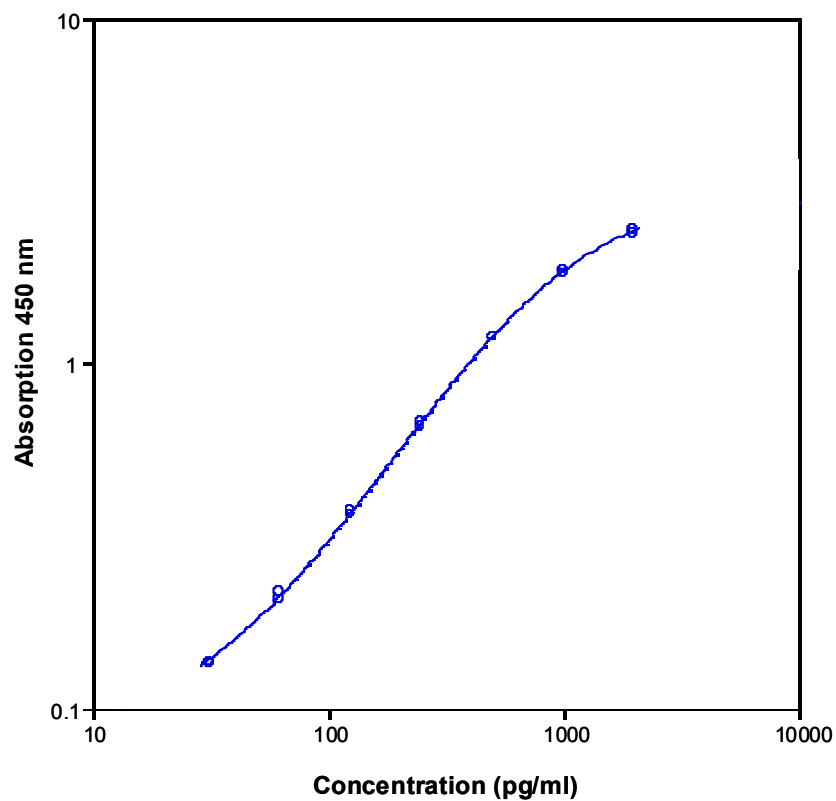

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

Table 2

Typical data using the pig IFN- γ ELISA

Measuring wavelength: 450 nm, Reference wavelength: 620 nm

Standard	Pig IFN- γ Concentration (pg/ml)	O.D. at 450 nm	Mean O.D. at 450 nm	C.V. (%)
1	2000	2.387	2.372	1.9
	2000	2.452		
2	1000	1.860	1.794	1.4
	1000	1.823		
3	500	1.198	1.145	0.7
	500	1.187		
4	250	0.655	0.620	2.6
	250	0.680		
5	125	0.377	0.324	2.1
	125	0.366		
6	63	0.217	0.164	3.3
	63	0.207		
7	31	0.135	0.088	0.5
	31	0.136		
Blank	0	0.050	0.048	
	0	0.046		

The OD values of the standard curve may vary according to the conditions of assay performance (e.g. operator, pipetting technique, washing technique or temperature effects). Furthermore shelf life of the kit may affect enzymatic activity and thus colour intensity. Values measured are still valid.

LIMITATIONS

- Since exact conditions may vary from assay to assay, a standard curve must be established for every run.
- Bacterial or fungal contamination of either screen samples or reagents or cross-contamination between reagents may cause erroneous results.
- Disposable pipette tips, flasks or glassware are preferred, reusable glassware must be washed and thoroughly rinsed of all detergents before use.
- Improper or insufficient washing at any stage of the procedure will result in either false positive or false negative results. Empty wells completely before dispensing fresh wash solution, fill with Wash Buffer as indicated for each wash cycle and do not allow wells to sit uncovered or dry for extended periods.

DRG® Interferon γ (porcine) (EIA-4828)**Revised 27 Jan. 2011 rm (Vers. 3.1)****For Veterinary Use Only****Sample Stability****1.1.2 Freeze-Thaw Stability**

Aliquots of spiked serum samples were stored at -20°C and thawed 5 times, and the pig IFN- γ levels determined. There was no significant loss of pig IFN- γ immunoreactivity detected by freezing and thawing.

1.1.3 Storage Stability

Aliquots of spiked serum samples were stored at -20°C, 2-8°C, room temperature (RT) and at 37°C, and the pig IFN- γ level determined after 24 h. There was no significant loss of pig IFN- γ immunoreactivity detected during storage under above conditions.

BIBLIOGRAPHY

1. Ciampolillo, A., Guastamacchia, E., Caragiulo, L., Lollino, G., De Robertis, O., Lattanzi, V., and Giorgino, R. (1993). In vitro secretion of interleukin-1 beta and interferon-gamma by peripheral blood lymphomononuclear cells in diabetic patients. *Diabetes Res. Clin. Pract.* 21, 87-93.
2. Edwards, B. S., Merritt, J. A., Fuhlbrigge, R. C., and Borden, E. C. (1985). Low doses of interferon alpha result in more effective clinical natural killer cell activation. *J. Clin. Invest.* 75, 1908-1913.
3. Gajewski, T. F., and Fitch, F. W. (1993). Anti-proliferative effect of IFN-gamma in immune regulation. I. IFN-gamma inhibits the proliferation of Th2 but not Th1 murine helper T lymphocyte clones. *J. Immunol.* 140, 4245-4252.
4. Le thi Bich Thuy, Queen, C., and Fauci, A. S. (1986). Interferon-gamma induces light chain synthesis in interleukin 2 stimulated human B cells. *Eur. J. Immunol.* 16, 547-550.
5. Mosmann, T. R., Cherwinski, H., Bond, M. W., Giedlin, M. A., and Coffman, R. L. (1986). Two types of murine helper T cell clone. I. Definition according to profiles of lymphokine activities and secreted proteins. *J. Immunol.* 136, 2348-2357.
6. Romagnani, S., Giudizi, M. G., Biagiotti, R., Almerigogna, F., Mingari, C., Maggi, E., Liang, C. M., and Moretta, L. (1986). B cell growth factor activity of interferon-gamma. Recombinant human interferon-gamma promotes proliferation of anti-mu-activated human B lymphocytes. *J. Immunol.* 136, 3513-3516.
7. Sastre, L., Roman, J. M., Teplow, D. B., Dreyer, W. J., Gee, C. E., Larson, R. S., Roberts, T. M., and Springer, T. A. (1986). A partial genomic DNA clone for the alpha subunit of the mouse complement receptor type 3 and cellular adhesion molecule Mac-1. *Proc. Natl. Acad. Sci. U. S. A.* 83, 5644-5648.
8. Sidman, C. L., Marshall, J. D., Shultz, L. D., Gray, P. W., and Johnson, H. M. (1984). Gamma-interferon is one of several direct B cell-maturing lymphokines. *Nature* 309, 801-804.
9. Urban, J. L., Shepard, H. M., Rothstein, J. L., Sugarman, B. J., and Schreiber, H. (1986). Tumor necrosis factor: a potent effector molecule for tumor cell killing by activated macrophages. *Proc. Natl. Acad. Sci. U. S. A.* 83, 5233-5237.

DRG® Interferon γ (porcine) (EIA-4828)

Revised 27 Jan. 2011 rm (Vers. 3.1)

For Veterinary Use Only

REAGENT PREPARATION SUMMARY

Wash Buffer (1x)

Add **Wash Buffer Concentrate** 20x (50 ml) to 950 ml distilled water.

Number of Strips	Wash Buffer Concentrate (ml)	Distilled Water (ml)
1 - 6	25	475
1 - 12	50	950

Assay Buffer (1x)

Add **Assay Buffer Concentrate** 20x (5 ml) to 95 ml distilled water.

Number of Strips	Assay Buffer Concentrate (ml)	Distilled Water (ml)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

Biotin-Conjugate

Make a 1:100 dilution of **Biotin-Conjugate** in Assay Buffer (1x):

Number of Strips	Biotin-Conjugate (ml)	Assay Buffer (1x) (ml)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

Streptavidin-HRP

Make a 1:100 dilution of **Streptavidin-HRP** in Assay Buffer (1x):

Number of Strips	Streptavidin-HRP (ml)	Assay Buffer (1x) (ml)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

Pig IFN- γ Standard

Reconstitute lyophilized **pig IFN- γ standard** with distilled water. (Reconstitution volume is stated on the label of the standard vial.)

DRG[®] Interferon γ (porcine) (EIA-4828)**Revised 27 Jan. 2011 rm (Vers. 3.1)****For Veterinary Use Only****TEST PROTOCOL SUMMARY**

1. Determine the number of microwell strips required.
2. Wash microwell strips twice with Wash Buffer.
3. Standard dilution on the microwell plate: Add 100 μ l Assay Buffer (1x), in duplicate, to all standard wells. Pipette 100 μ l prepared standard into the first wells and create standard dilutions by transferring 100 μ l from well to well. Discard 100 μ l from the last wells.
Alternatively external standard dilution in tubes (see 1.1.1): Pipette 100 μ l of these standard dilutions in the microwell strips.
4. Add 100 μ l Assay Buffer (1x), in duplicate, to the blank wells.
5. Add 50 μ l Assay Buffer (1x) to sample wells.
6. Add 50 μ l sample in duplicate, to designated sample wells.
7. Cover microwell strips and incubate 1 hour at room temperature (18° to 25°C).
8. Prepare Biotin-Conjugate.
9. Empty and wash microwell strips 3 times with Wash Buffer.
10. Add 100 μ l Biotin-Conjugate to all wells.
11. Cover microwell strips and incubate 1 hour at room temperature (18° to 25°C).
12. Prepare Streptavidin-HRP.
13. Empty and wash microwell strips 3 times with Wash Buffer.
14. Add 100 μ l diluted Streptavidin-HRP to all wells.
15. Cover microwell strips and incubate 30 minutes at room temperature (18° to 25°C).
16. Empty and wash microwell strips 3 times with Wash Buffer.
17. Add 100 μ l of TMB Substrate Solution to all wells.
18. Incubate the microwell strips for about 10 minutes at room temperature (18° to 25°C).
19. Add 100 μ l Stop Solution to all wells.
20. Blank microwell reader and measure colour intensity at 450 nm.

Note: If instructions in this protocol have been followed samples have been diluted 1:2 (50 μ l sample + 50 μ l Assay Buffer (1x)), the concentration read from the standard curve must be multiplied by the dilution factor (x 2).