

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

INTENDED USE

The mouse Interleukin-12 (Mouse) ELISA is an enzyme-linked immunosorbent assay for the quantitative detection of mouse Interleukin-12 (IL-12).

The mouse IL-12 ELISA is for research use only. Not for diagnostic or therapeutic procedures.

SUMMARY

Interleukin-12 (IL-12) is a pleiotropic cytokine, formerly termed cytotoxic lymphocyte maturation factor (CLMF) or natural killer cell stimulatory factor (NKSF) (1,3,4), which is produced primarily by stimulated macrophages. IL-12 has been shown to be a proinflammatory cytokine produced by phagocytic cells (5), B cells (1,3), and other antigen - presenting cells that modulate adaptive immune responses by favoring the generation of T-helper type 1 cells (2).

IL-12 exerts a variety of biological effects on T and natural killer cells. Apart from promotion of Th1 development and its ability to promote cytolytic activity it mediates some of its physiological activities by acting as a potent inducer of interferon gamma (IFN γ) production and the stimulation of other cytokines from peripheral blood T and NK cells, (6, 7). IFN γ then enhances the ability of the phagocytic cells to produce IL-12 and other proinflammatory cytokines. Thus, IL-12 induced IFN γ acts in a positive feedback loop that represents an important amplifying mechanism in the inflammatory response to infections (2).

Its role in directing development of a Th1 type immune response from naive T cells demonstrates its critical role in regulation of the immune response and strongly suggests its potential usefulness in cancer therapy (4).

PRINCIPLES OF THE TEST

An anti-mouse IL-12 coating antibody is adsorbed onto microwells.

Mouse IL-12 present in the sample or standard binds to antibodies adsorbed to the microwells. A biotin-conjugated anti-mouse IL-12 antibody is added and binds to mouse IL-12 captured by the first antibody.

Following incubation unbound biotin-conjugated anti-mouse IL-12 antibody is removed during a wash step. Streptavidin-HRP is added and binds to the biotin-conjugated anti-mouse IL-12 antibody.

Following incubation unbound Streptavidin-HRP is removed during a wash step, and substrate solution reactive with HRP is added to the wells.

A coloured product is formed in proportion to the amount of mouse IL-12 present in the sample or standard. The reaction is terminated by addition of acid and absorbance is measured at 450 nm. A standard curve is prepared from 7 mouse IL-12 standard dilutions and mouse IL-12 sample concentration determined.

REAGENTS PROVIDED

Reagents for mouse IL-12 ELISA (96 tests)

- 1 aluminium pouch with a **Microwell Plate coated** with monoclonal antibody to mouse IL-12
- 1 vial (100 μ L) **Biotin-Conjugate** anti-mouse IL-12 monoclonal antibody
- 1 vial (150 μ L) **Streptavidin-HRP**

DRG International Inc., USA

Fax: (908) 233-0758 • E-mail: corp@drg-international.com • Web: www.drg-international.com

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

- 2 vials mouse IL-12 **Standard** lyophilized, 4000 pg/mL upon reconstitution
- 1 vial (5 mL) **Assay Buffer Concentrate** 20x (PBS with 1% Tween 20 and 10% BSA)
- 1 bottle (50 mL) **Wash Buffer Concentrate** 20x (PBS with 1% Tween 20)
- 1 vial (7 mL) **Substrate Solution I** (tetramethyl-benzidine)
- 1 vial (7 mL) **Substrate Solution II** (0.02 % buffered hydrogen peroxide)
- 1 vial (12 mL) **Stop Solution** (1M Phosphoric acid)
- 1 vial (0.4 mL) **Blue-Dye**
- 1 vial (0.4 mL) **Green-Dye**
- 1 vial (0.4 mL) **Red-Dye**
- 4 **Adhesive Films**

STORAGE INSTRUCTIONS – ELISA KIT

Store kit reagents between 2°C and 8°C. Immediately after use remaining reagents should be returned to cold storage (2°C to 8°C). Expiry of the kit and reagents is stated on labels.

Expiry of the kit components can only be guaranteed if the components are stored properly, and if, in case of repeated use of one component, this reagent is not contaminated by the first handling.

SPECIMEN COLLECTION AND STORAGE INSTRUCTIONS

Cell culture supernatant and serum were tested with this assay. Other body fluids might be suitable for use in the assay. Remove serum from the clot as soon as possible after clotting.

Pay attention to a possible “**Hook Effect**” due to high sample concentrations (see 0).

Samples containing a visible precipitate must be clarified prior to use in the assay. Do not use grossly hemolyzed or lipemic specimens.

Samples should be aliquoted and must be stored frozen at -20°C to avoid loss of bioactive mouse IL-12. If samples are to be run within 24 hours, they may be stored at 2°C to 8°C (for sample stability refer to 1.12).

Avoid repeated freeze-thaw cycles. Prior to assay, the frozen sample should be brought to room temperature slowly and mixed gently.

MATERIALS REQUIRED BUT NOT PROVIDED

- 5 mL and 10 mL graduated pipettes
- 5 µL to 1000 µL adjustable single channel micropipettes with disposable tips
- 50 µL to 300 µL adjustable multichannel micropipette with disposable tips

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

- Multichannel micropipette reservoir
- Beakers, flasks, cylinders necessary for preparation of reagents
- Device for delivery of wash solution (multichannel wash bottle or automatic wash system)
- Microwell strip reader capable of reading at 450 nm (620 nm as optional reference wave length)
- Glass-distilled or deionized water
- Statistical calculator with program to perform regression analysis

PRECAUTIONS FOR USE

- All chemicals should be considered as potentially hazardous. We therefore recommend that this product is handled only by those persons who have been trained in laboratory techniques and that it is used in accordance with the principles of good laboratory practice. Wear suitable protective clothing such as laboratory overalls, safety glasses and gloves. Care should be taken to avoid contact with skin or eyes. In the case of contact with skin or eyes wash immediately with water. See material safety data sheet(s) and/or safety statement(s) for specific advice.
- Reagents are intended for research use only and are not for use in diagnostic or therapeutic procedures.
- Do not mix or substitute reagents with those from other lots or other sources.
- Do not use kit reagents beyond expiration date on label.
- Do not expose kit reagents to strong light during storage or incubation.
- Do not pipette by mouth.
- Do not eat or smoke in areas where kit reagents or samples are handled.
- Avoid contact of skin or mucous membranes with kit reagents or specimens.
- Rubber or disposable latex gloves should be worn while handling kit reagents or specimens.
- Avoid contact of substrate solutions with oxidizing agents and metal.
- Avoid splashing or generation of aerosols.
- In order to avoid microbial contamination or cross-contamination of reagents or specimens which may invalidate the test use disposable pipette tips and/or pipettes.
- Use clean, dedicated reagent trays for dispensing the conjugate and substrate reagent.
- Exposure to acid inactivates the conjugate.
- Glass-distilled water or deionized water must be used for reagent preparation.
- Substrate solutions must be at room temperature prior to use.
- Decontaminate and dispose specimens and all potentially contaminated materials as they could contain infectious agents. The preferred method of decontamination is autoclaving for a minimum of 1 hour at 121.5°C.
- Liquid wastes not containing acid and neutralized waste may be mixed with sodium hypochlorite in volumes such that the final mixture contains 1.0% sodium hypochlorite. Allow 30 minutes for effective decontamination. Liquid waste containing acid must be neutralized prior to the addition of sodium hypochlorite.

PREPARATION OF REAGENTS

Buffer concentrates should be brought to room temperature and should be diluted before starting the test procedure. If crystals have formed in **buffer concentrates**, warm them gently until crystals have completely dissolved.

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

1.1 Wash Buffer (1x)

If crystals have formed in the **Wash Buffer Concentrate** (20x), warm it gently until they have completely dissolved. Pour entire contents (50 mL) of the Wash Buffer Concentrate into a clean 1000 mL graduated cylinder. Bring to final volume of 1000 mL with glass-distilled or deionized water. Mix gently to avoid foaming. The pH of the final solution should adjust to 7.4.

Transfer to a clean wash bottle and store at 2°C to 25°C. Please note that Wash Buffer (1x) is stable for 30 days.

Wash Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Wash Buffer Concentrate (mL)	Distilled Water (mL)
1 - 6	25	475
1 - 12	50	950

1.2 Assay Buffer (1x)

Pour the entire contents (5 mL) of the **Assay Buffer Concentrate** (20x) into a clean 100 mL graduated cylinder. Bring to final volume of 100 mL with distilled water. Mix gently to avoid foaming.

Store at 2°C to 8°C. Please note that the Assay Buffer (1x) is stable for 30 days.

Assay Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Assay Buffer Concentrate (mL)	Distilled Water (mL)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

1.3 Biotin-Conjugate

Please note that the Biotin-Conjugate should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated **Biotin-Conjugate** solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Biotin-Conjugate (mL)	Assay Buffer (1x) (mL)
1 - 6	0.03	2.97
1 - 12	0.06	5.94

1.4 Streptavidin-HRP

Please note that the Streptavidin-HRP should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated **Streptavidin-HRP** solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

DRG International Inc., USA

Fax: (908) 233-0758 • E-mail: corp@drg-international.com • Web: www.drg-international.com

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

Number of Strips	Streptavidin-HRP (mL)	Assay Buffer (1x) (mL)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

1.5 Mouse IL-12 Standard

Reconstitute **mouse IL-12 standard** by addition of distilled water.

Reconstitution volume is stated on the label of the standard vial. Swirl or mix gently to insure complete and homogeneous solubilisation (concentration of reconstituted standard = 4000 pg/mL).

The standard has to be used immediately after reconstitution and cannot be stored.

Standard dilutions can be prepared directly on the microwell plate (see 10.c) or alternatively in tubes (see 1.5.1).

1.5.1 External Standard Dilution

Label 7 tubes, one for each standard point: S1, S2, S3, S4, S5, S6, S7

Then prepare 1:2 serial dilutions for the standard curve as follows:

Pipette 225 µL of Assay Buffer (1x) into each tube.


Pipette 225 µL of reconstituted standard (concentration = 4000 pg/mL) into the first tube, labelled S1, and mix (concentration of standard 1 = 2000 pg/mL).

Pipette 225 µL of this dilution into the second tube, labelled S2, and mix thoroughly before the next transfer.

Repeat serial dilutions 5 more times thus creating the points of the standard curve (see Figure 1).

Assay Buffer (1x) serves as blank.

Figure 1

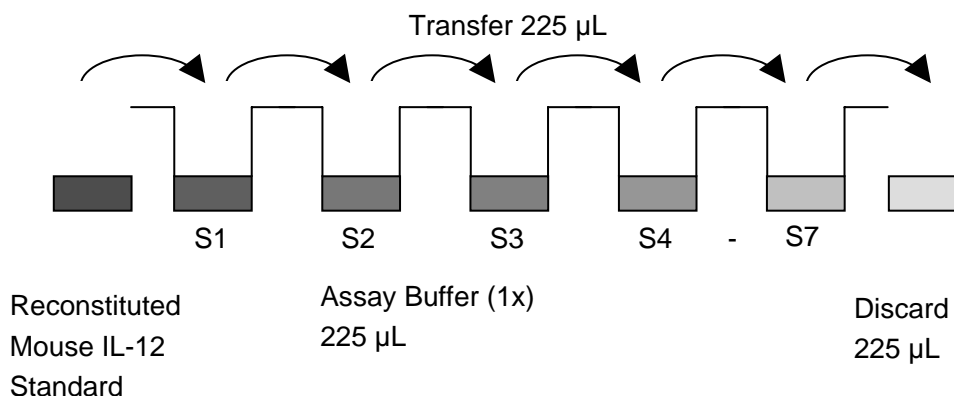

DRG® Interleukin 12 (Mouse) (EIA-4920)**As of 14 Nov. 2008 (Vers. 1.0)****For Veterinary Use Only****1.6 TMB Substrate Solution**

Using clean pipettes and containers known to be metal free, dispense an equal volume of **Substrate Solution I** into **Substrate Solution II** and swirl gently to mix.

The TMB Substrate Solution may develop a yellow tinge over time. This does not seem to affect product performance. A blue colour present in the TMB Substrate Solution, however, indicates that it has been contaminated and must be discarded.

The TMB Substrate Solution must be used within a few minutes after mixing.

Warm to room temperature before use.

Avoid direct exposure of TMB reagents to intense light and oxidizing agents during storage or incubation.

Substrate preparation according to assay size:

Number of Strips	Substrate Solution I (mL)	Substrate Solution II (mL)
1 - 6	3.0	3.0
1 - 12	6.0	6.0

1.7 Addition of Colour-giving Reagents: Blue-Dye, Green-Dye, Red-Dye

In order to help our customers to avoid any mistakes in pipetting the ELISAs, a tool is offered that helps to monitor the addition of even very small volumes of a solution to the reaction well by giving distinctive colours to each step of the ELISA procedure.

This procedure is optional, does not in any way interfere with the test results, and is designed to help the customer with the performance of the test, but can also be omitted, just following the instruction booklet.

Alternatively, the dye solutions from the stocks provided (**Blue-Dye, Green-Dye, Red-Dye**) can be added to the reagents according to the following guidelines:

1. Diluent: Before standard and sample dilution add the **Blue-Dye** at a dilution of 1:250 (see table below) to the appropriate diluent (1x) according to the test protocol. After addition of **Blue-Dye**, proceed according to the instruction booklet.

5 mL Assay Buffer (1x)	20 µL Blue-Dye
12 mL Assay Buffer (1x)	48 µL Blue-Dye
50 mL Assay Buffer (1x)	200 µL Blue-Dye

2. Biotin-Conjugate: Before dilution of the concentrated Biotin-Conjugate, add the **Green-Dye** at a dilution of 1:100 (see table below) to the Assay Buffer (1x) used for the final conjugate dilution. Proceed after addition of **Green-Dye** according to the instruction booklet: Preparation of Biotin-Conjugate.

3 mL Assay Buffer (1x)	30 µL Green-Dye
6 mL Assay Buffer (1x)	60 µL Green-Dye

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only


3. Streptavidin-HRP: Before dilution of the concentrated Streptavidin-HRP, add the **Red-Dye** at a dilution of 1:250 (see table below) to the Assay Buffer (1x) used for the final Streptavidin-HRP dilution. Proceed after addition of **Red-Dye** according to the instruction booklet: Preparation of Streptavidin-HRP.

6 mL Assay Buffer (1x)	24 µL Red-Dye
12 mL Assay Buffer (1x)	48 µL Red-Dye

TEST PROTOCOL

- Determine the number of microwell strips required to test the desired number of samples plus appropriate number of wells needed for running blanks and standards. Each sample, standard, blank and optional control sample should be assayed in duplicate. Remove extra microwell strips from holder and store in foil bag with the desiccant provided at 2°C-8°C sealed tightly.
- Wash the microwell strips twice with approximately 400 µL **Wash Buffer** per well with thorough aspiration of microwell contents between washes. Allow the Wash Buffer to sit in the wells for about **10 – 15 seconds** before aspiration. Take care not to scratch the surface of the microwells.
After the last wash step, empty wells and tap microwell strips on absorbent pad or paper towel to remove excess Wash Buffer. Use the microwell strips immediately after washing. Alternatively microwell strips can be placed upside down on a wet absorbent paper for not longer than 15 minutes. **Do not allow wells to dry.**
- Standard dilution on the microwell plate** (Alternatively the standard dilution can be prepared in tubes - see 1.5.1): Add 100 µL of Assay Buffer (1x) in duplicate to all **standard wells**. Pipette 100 µL of prepared **standard** (see Preparation of Standard 1.5, concentration = 4000 pg/mL) in duplicate into well A1 and A2 (see Table 1). Mix the contents of wells A1 and A2 by repeated aspiration and ejection (concentration of standard 1, S1 = 2000 pg/mL), and transfer 100 µL to wells B1 and B2, respectively (see Figure 2). Take care not to scratch the inner surface of the microwells. Continue this procedure 5 times; creating two rows of mouse IL-12 standard dilutions ranging from 2000.0 to 31.3 pg/mL. Discard 100 µL of the contents from the last microwells (G1, G2) used.

Figure 2

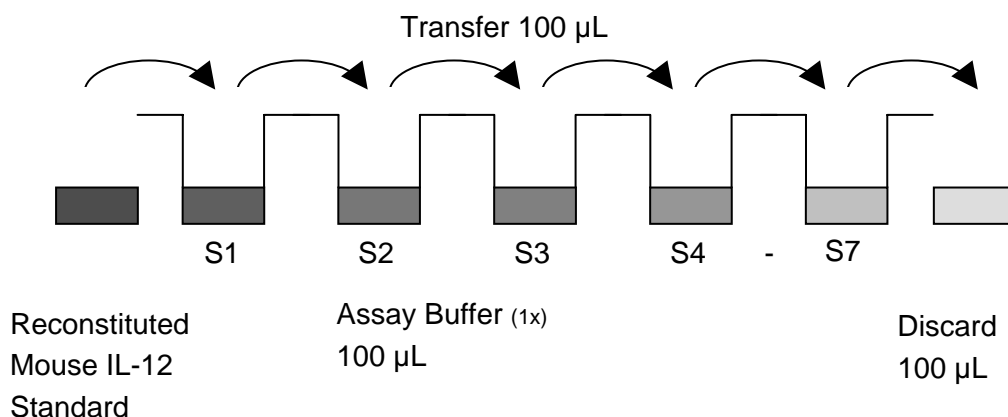

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

In case of an **external standard dilution** (see 1.5.1), pipette 100 µL of these standard dilutions (S1 - S7) in the standard wells according to Table 1.

Table 1:

Table depicting an example of the arrangement of blanks, standards and samples in the microwell strips:

	1	2	3	4
A	Standard 1 (2000.0 pg/mL)	Standard 1 (2000.0 pg/mL)	Sample 1	Sample 1
B	Standard 2 (1000.0 pg/mL)	Standard 2 (1000.0 pg/mL)	Sample 2	Sample 2
C	Standard 3 (500.0 pg/mL)	Standard 3 (500.0 pg/mL)	Sample 3	Sample 3
D	Standard 4 (250.0 pg/mL)	Standard 4 (250.0 pg/mL)	Sample 4	Sample 4
E	Standard 5 (125.0 pg/mL)	Standard 5 (125.0 pg/mL)	Sample 5	Sample 5
F	Standard 6 (62.5 pg/mL)	Standard 6 (62.5 pg/mL)	Sample 6	Sample 6
G	Standard 7 (31.3 pg/mL)	Standard 7 (31.3 pg/mL)	Sample 7	Sample 7
H	Blank	Blank	Sample 8	Sample 8

- d. Add 100 µL of **Assay Buffer (1x)** in duplicate to the **blank wells**.
- e. Add 50 µL of **Assay Buffer (1x)** to the **sample wells**.
- f. Add 50 µL of each **sample** in duplicate to the **sample wells**.
- g. Prepare **Biotin-Conjugate** (see Preparation of Biotin-Conjugate 1.3).
- h. Add 50 µL of **Biotin-Conjugate** to all wells.
- i. Cover with an adhesive film and incubate at room temperature (18°C to 25°C) for 2 hours, if available on a microplate shaker set at 200 rpm.
- j. Prepare **Streptavidin-HRP** (refer to Preparation of Streptavidin-HRP 0).
- k. Remove adhesive film and empty wells. **Wash** microwell strips 3 times according to point b. of the test protocol. Proceed immediately to the next step.
- l. Add 100 µL of diluted **Streptavidin-HRP** to all wells, including the blank wells.
- m. Cover with an adhesive film and incubate at room temperature (18°C to 25°C) for 1 hour, if available on a microplate shaker set at 200 rpm.
- n. Prepare **TMB Substrate Solution** a few minutes prior to use. (refer to Preparation of TMB Substrate Solution 1.6).

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

- o. Remove adhesive film and empty wells. **Wash** microwell strips 3 times according to point b. of the test protocol. Proceed immediately to the next step.
- p. Pipette 100 µL of mixed **TMB Substrate Solution** to all wells.
- q. Incubate the microwell strips at room temperature (18°C to 25°C) for about 10 min. Avoid direct exposure to intense light.

The colour development on the plate should be monitored and the substrate reaction stopped (see next point of this protocol) before positive wells are no longer properly recordable.

Determination of the ideal time period for colour development has to be done individually for each assay.

It is recommended to add the stop solution when the highest standard has developed a dark blue colour. Alternatively the colour development can be monitored by the ELISA reader at 620 nm. The substrate reaction should be stopped as soon as Standard 1 has reached an OD of 0.6 – 0.65.

- r. Stop the enzyme reaction by quickly pipetting 100 µL of **Stop Solution** into each well. It is important that the Stop Solution is spread quickly and uniformly throughout the microwells to completely inactivate the enzyme. Results must be read immediately after the Stop Solution is added or within one hour if the microwell strips are stored at 2 - 8°C in the dark.
- s. Read absorbance of each microwell on a spectro-photometer using 450 nm as the primary wave length (optionally 620 nm as the reference wave length; 610 nm to 650 nm is acceptable). Blank the plate reader according to the manufacturer's instructions by using the blank wells. Determine the absorbance of both the samples and the standards.

Note: In case of incubation without shaking the obtained O.D. values may be lower than indicated below. Nevertheless the results are still valid.

CALCULATION OF RESULTS

- Calculate the average absorbance values for each set of duplicate standards and samples. Duplicates should be within 20 per cent of the mean value.
- Create a standard curve by plotting the mean absorbance for each standard concentration on the ordinate against the mouse IL-12 concentration on the abscissa. Draw a best fit curve through the points of the graph (a 5-parameter curve fit is recommended).
- To determine the concentration of circulating mouse IL-12 for each sample, first find the mean absorbance value on the ordinate and extend a horizontal line to the standard curve. At the point of intersection, extend a vertical line to the abscissa and read the corresponding mouse IL-12 concentration.
- **If instructions in this protocol have been followed samples have been diluted 1:2 (50 µL sample + 50 µL Assay Buffer (1x)), the concentration read from the standard curve must be multiplied by the dilution factor (x 2).**
- **Calculation of samples with a concentration exceeding standard 1 may result in incorrect, low mouse IL-12 levels (Hook Effect). Such samples require further external predilution according to expected mouse IL-12 values with Assay Buffer (1x) in order to precisely quantitate the actual mouse IL-12 level.**
- It is suggested that each testing facility establishes a control sample of known mouse IL-12 concentration and runs this additional control with each assay. If the values obtained are not within the expected range of the control, the assay results may be invalid.

DRG® Interleukin 12 (Mouse) (EIA-4920)


As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

- A representative standard curve is shown in Figure 3. This curve cannot be used to derive test results. Each laboratory must prepare a standard curve for each group of microwell strips assayed.

Figure 3:

Representative standard curve for mouse IL-12 ELISA. Mouse IL-12 was diluted in serial 2-fold steps in Assay Buffer (1x). Do not use this standard curve to derive test results. A standard curve must be run for each group of microwell strips assayed.

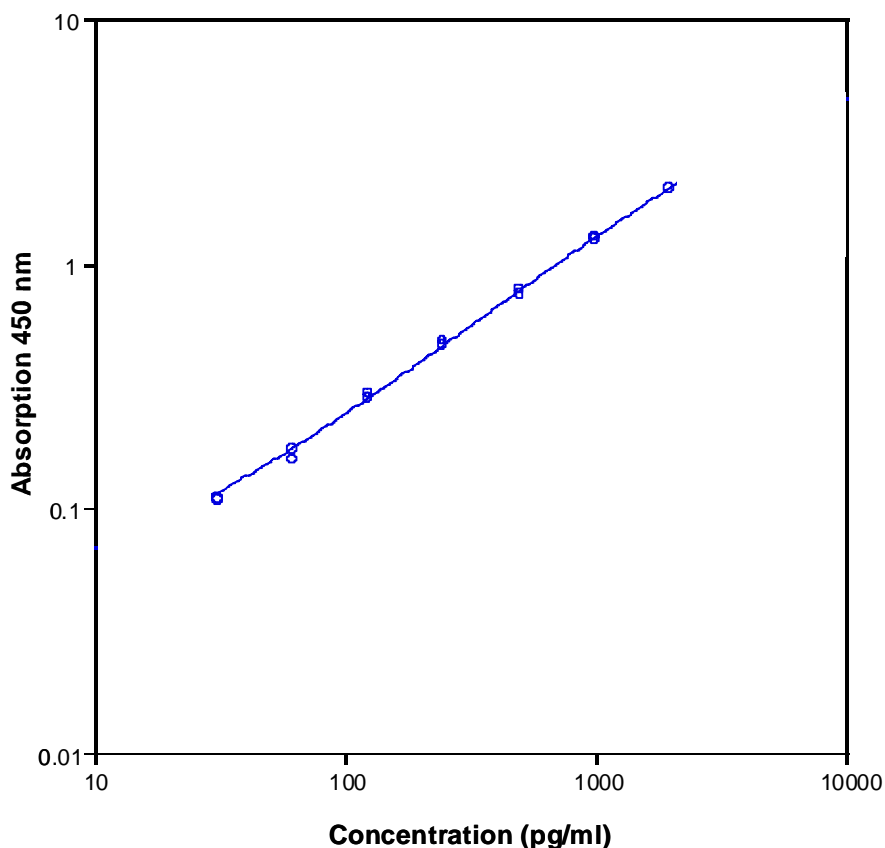

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

Table 2: Typical data using the mouse IL-12 ELISA
Measuring wavelength: 450 nm, Reference wavelength: 620 nm

Standard	Mouse IL-12 Concentration (pg/mL)	O.D. at 450 nm	Mean O.D. at 450 nm	C.V. (%)
1	2000.0	2.029	2.019	0.2
	2000.0	2.020		
2	1000.0	1.242	1.259	1.8
	1000.0	1.287		
3	500.0	0.787	0.762	2.5
	500.0	0.748		
4	250.0	0.462	0.466	1.9
	250.0	0.480		
5	125.0	0.281	0.281	1.9
	125.0	0.292		
6	62.5	0.158	0.160	4.5
	62.5	0.173		
7	31.3	0.108	0.109	0.5
	31.3	0.109		
Blank	0	0.041	0.046	
	0	0.050		

The OD values of the standard curve may vary according to the conditions of assay performance (e.g. operator, pipetting technique, washing technique or temperature effects). Furthermore shelf life of the kit may affect enzymatic activity and thus colour intensity. Values measured are still valid.

LIMITATIONS

- Since exact conditions may vary from assay to assay, a standard curve must be established for every run.
- Bacterial or fungal contamination of either screen samples or reagents or cross-contamination between reagents may cause erroneous results.
- Disposable pipette tips, flasks or glassware are preferred, reusable glassware must be washed and thoroughly rinsed of all detergents before use.
- Improper or insufficient washing at any stage of the procedure will result in either false positive or false negative results. Empty wells completely before dispensing fresh wash solution, fill with Wash Buffer as indicated for each wash cycle and do not allow wells to sit uncovered or dry for extended periods.

DRG® Interleukin 12 (Mouse) (EIA-4920)**As of 14 Nov. 2008 (Vers. 1.0)****For Veterinary Use Only****PERFORMANCE CHARACTERISTICS****1.8 Sensitivity**

The limit of detection of mouse IL-12 defined as the analyte concentration resulting in an absorbance significantly higher than that of the dilution medium (mean plus 2 standard deviations) was determined to be 4.0 pg/mL (mean of 6 independent assays).

1.9 Reproducibility**1.9.1 Intra-assay**

Reproducibility within the assay was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 8 serum samples containing different concentrations of mouse IL-12. 2 standard curves were run on each plate. The calculated overall intra-assay coefficient of variation was < 5%.

1.9.2 Inter-assay

Assay to assay reproducibility within one laboratory was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 8 serum samples containing different concentrations of mouse IL-12. 2 standard curves were run on each plate. The calculated overall inter-assay coefficient of variation was < 10%.

1.10 Spike Recovery

The spike recovery was evaluated by spiking 4 levels of mouse IL-12 into pooled murine serum. Recoveries were determined in 2 independent experiments with 4 replicates each.

The unspiked serum was used as blank in these experiments.

The overall mean recovery was 93%.

1.11 Dilution Parallelism

4 serum samples with different levels of mouse IL-12 were analysed at serial 2 fold dilutions with 4 replicates each.

The overall mean recovery was 104%.

1.12 Sample Stability**1.12.1 Freeze-Thaw Stability**

Aliquots of spiked serum samples were stored at -20°C and thawed 5 times, and the mouse IL-12 levels determined. There was no significant loss of mouse IL-12 immunoreactivity detected by freezing and thawing.

1.12.2 Storage Stability

Aliquots of spiked serum samples were stored at -20°C, 2-8°C, room temperature (RT) and at 37°C, and the mouse IL-12 level determined after 24 h. There was no significant loss of mouse IL-12 immunoreactivity detected during storage at -20°C, 2-8°C and RT.

A significant loss of mouse IL-12 immunoreactivity (50%) was detected during storage at 37°C after 24 h.

DRG® Interleukin 12 (Mouse) (EIA-4920)**As of 14 Nov. 2008 (Vers. 1.0)****For Veterinary Use Only****1.13 Specificity**

The interference of circulating factors of the immune system was evaluated by spiking these proteins at physiologically relevant concentrations into mouse IL-12 positive serum.

There was no cross reactivity detected.

1.14 Expected Values

There were no detectable mouse IL-12 levels found.

Elevated mouse IL-12 levels depend on the type of immunological disorder.

BIBLIOGRAPHY

1. Kobayashi et al., (1989). J. Exp. Med., 170:827.
2. Ma X., Chow J.M., Gri G., Carra G., Gerosa F., Wolf S.F., Dzialo R., Trinchieri G., (1996). The interleukin-12 p40 gene promoter is primed by interferon gamma in monocytic cells. J. Exp. Med. 183(1):147-57.
3. Stern et al., (1990). Proc. Natl. Acad. Sci. 87:6808.
4. Tahara H., Lotze M.T., (1995). Antitumor effects of interleukin-12 (IL-12): applications for the immunotherapy and gene therapy of cancer. Gene Ther. 2(2):96-106.
5. Trinchieri et al., (1992). Prog. Growth Factor Res., 4:355.
6. Wolf S.F., Sieburth D., Sypek J., (1994). Interleukin-12: a key modulator of immune function. Stem Cells Dayt. 12(2):154-68.
7. Zeh, H.J. et al (1994): The Cytokine Handbook, 2nd edition, Thomson, A. edition, Academic Press, New York, P239.

REAGENT PREPARATION SUMMARY**1.15 Wash Buffer (1x)**

Add **Wash Buffer Concentrate** 20x (50 mL) to 950 mL distilled water.

Number of Strips	Wash Buffer Concentrate (mL)	Distilled Water (mL)
1 - 6	25	475
1 - 12	50	950

1.16 Assay Buffer (1x)

Add **Assay Buffer Concentrate** 20x (5 mL) to 95 mL distilled water.

Number of Strips	Assay Buffer Concentrate (mL)	Distilled Water (mL)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

1.17 Biotin-Conjugate

Make a 1:100 dilution of **Biotin-Conjugate** in Assay Buffer (1x):

DRG® Interleukin 12 (Mouse) (EIA-4920)**As of 14 Nov. 2008 (Vers. 1.0)****For Veterinary Use Only**

Number of Strips	Biotin-Conjugate (mL)	Assay Buffer (1x) (mL)
1 - 6	0.03	2.97
1 - 12	0.06	5.94

1.18 Streptavidin-HRPMake a 1:100 dilution of **Streptavidin-HRP** in Assay Buffer (1x):

Number of Strips	Streptavidin-HRP (mL)	Assay Buffer (1x) (mL)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

1.19 Mouse IL-12 StandardReconstitute lyophilized **mouse IL-12 standard** with distilled water. (Reconstitution volume is stated on the label of the standard vial.)**1.20 TMB Substrate Solution**Dispense an equal volume of **Substrate Solution I** into **Substrate Solution II**.

Number of Strips	Substrate Solution I (mL)	Substrate Solution II (mL)
1 - 6	3.0	3.0
1 - 12	6.0	6.0

TEST PROTOCOL SUMMARY

1. Determine the number of microwell strips required.
2. Wash microwell strips twice with Wash Buffer.
3. Standard dilution on the microwell plate: Add 100 µL Assay Buffer (1x), in duplicate, to all standard wells. Pipette 100 µL prepared standard into the first wells and create standard dilutions by transferring 100 µL from well to well. Discard 100 µL from the last wells.
Alternatively external standard dilution in tubes (see 1.5.1): Pipette 100 µL of these standard dilutions in the microwell strips.
4. Add 100 µL Assay Buffer (1x), in duplicate, to the blank wells.
5. Add 50 µL Assay Buffer (1x) to sample wells.
6. Add 50 µL sample in duplicate, to designated sample wells.
7. Prepare Biotin-Conjugate.
8. Add 50 µL Biotin-Conjugate to all wells.
9. Cover microwell strips and incubate 2 hours at room temperature (18° to 25°C).

DRG[®] Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

For Veterinary Use Only

10. Prepare Streptavidin-HRP.
11. Empty and wash microwell strips 3 times with Wash Buffer.
12. Add 100 µL diluted Streptavidin-HRP to all wells.
13. Cover microwell strips and incubate 1 hour at room temperature (18° to 25°C).
14. Prepare TMB Substrate Solution few minutes prior to use.
15. Empty and wash microwell strips 3 times with Wash Buffer.
16. Add 100 µL of mixed TMB Substrate Solution to all wells.
17. Incubate the microwell strips for about 10 minutes at room temperature (18°to 25°C).
18. Add 100 µL Stop Solution to all wells.
19. Blank microwell reader and measure colour intensity at 450 nm.





















Note: If instructions in this protocol have been followed samples have been diluted 1:2 (50 µL sample + 50 µL Assay Buffer (1x)), the concentration read from the standard curve must be multiplied by the dilution factor (x 2).

DRG® Interleukin 12 (Mouse) (EIA-4920)

As of 14 Nov. 2008 (Vers. 1.0)

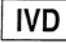For Veterinary Use Only

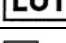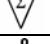SYMBOLS USED WITH DRG ASSAYS

Symbol	English	Deutsch	Français	Español	Italiano
	Consult instructions for use	Gebrauchsanweisung beachten	Consulter les instructions d'utilisation	Consulte las instrucciones de uso	Consultare le istruzioni per l'uso
	European Conformity	CE-Konformitätskennzeichnung	Conformité aux normes européennes	Conformidad europea	Conformità europea
	In vitro diagnostic device	In-vitro-Diagnostikum	Usage Diagnostic in vitro	Para uso Diagnóstico in vitro	Per uso Diagnostica in vitro
	For research use only	Nur für Forschungszwecke	Seulement dans le cadre de recherches	Sólo para uso en investigación	Solo a scopo di ricerca
	Catalogue number	Katalog-Nr.	Numéro de catalogue	Número de catálogo	Numero di Catalogo
	Lot. No. / Batch code	Chargen-Nr.	Numéro de lot	Número de lote	Numero di lotto
	Contains sufficient for <n> tests/	Ausreichend für "n" Ansätze	Contenu suffisant pour "n" tests	Contenido suficiente para <n> ensayos	Contenuto sufficiente per "n" saggi
	Storage Temperature	Lagerungstemperatur	Température de conservation	Temperatura de conservación	Temperatura di conservazione
	Expiration Date	Mindesthaltbarkeitsdatum	Date limite d'utilisation	Fecha de caducidad	Data di scadenza
	Legal Manufacturer	Hersteller	Fabricant	Fabricante	Fabbricante
Distributed by	Distributor	Vertreiber	Distributeur	Distribuidor	Distributore
Content	Content	Inhalt	Conditionnement	Contenido	Contenuto
Volume/No.	Volume / No.	Volumen/Anzahl	Volume/Quantité	Volumen/Número	Volume/Quantità
Symbol	Portugues	Dansk	Svenska	Ελληνικά	
	Consulte as instruções de utilização	Se brugsanvisning	Se bruksanvisningen	Εγχειρίδιο χρήστη	
	Conformidade com as normas europeias	Europeisk overensstemmelse	Europeisk överensstämmelse	Ευρωπαϊκή Συμμόρφωση	
	Diagnóstico in vitro	In vitro diagnostik	Diagnostik in vitro	in vitro διαγνωστικό	
					
	Catálogo n.º	Katalognummer	Katalog nummer	Αριθμός καταλόγου	
	No do lote	Lot nummer	Batch-nummer	Αριθμός Παρτίδος	
		Indeholder tilstrækkeligt til "n" test	Innehåller tillräckligt till "n" tester	Περιεχόμενο επαρκές για «n» εξετάσεις	
	Temperatura de conservação	Opbevarings-temperatur	Förvaringstemperatur	Θερμοκρασία αποθήκευσης	
	Prazo de validade	Udløbsdato	Bäst före datum	Ημερομηνία λήξης	
	Fabricante	Producent	Tillverkare	Κατασκευαστής	
Distributed by					
Content	Conteúdo	Indhold	Innehåll	Περιεχόμενο	
Volume/No.	Volume/Número	Volumen/antal	Volym/antal	Όγκος/αριθ..	