

RAT/MOUSE GROWTH HORMONE ELISA KIT 96-Well Plate (Cat. EZRMGH-45K)

I.	Intended Use	2
II.	Principles Of Procedure	2
III.	Reagents Supplied	2
IV.	Storage and Stability	4
۷.	Reagent Precautions	4
VI.	Materials Required But Not Provided	4
VII.	Sample Collection And Storage	5
VIII.	Sample Preparation	5
IX.	Standard and Quality Controls Preparation	6
Х.	Assay Procedure	7
XI.	Microtiter Plate Arrangement	10
XII.	Calculations	11
XIII.	Interpretation	11
XIV.	Standard Curve	12
XV.	Assay Characteristics	13
XVI.	Quality Controls	15
XVII.	Troubleshooting Guide	16
XVIII.	1 5	16
XIX.	Ordering Information	17

RAT/MOUSE GROWTH HORMONE ELISA KIT 96-Well Plate (Cat. #EZRMGH-45K)

I. INTENDED USE

This Rat/Mouse Growth Hormone ELISA kit is used for the non-radioactive quantification of Growth Hormone in rat or mouse serum, plasma, tissue extracts or cell culture media samples. One kit is sufficient to measure 39 unknown samples in duplicate. *This kit is for research purpose only.*

II. PRINCIPLES OF PROCEDURE

This assay is a Sandwich ELISA based, sequentially, on: 1) capture of rat or mouse Growth Hormone molecules from samples to the wells of a microtiter plate coated by a pre-titered amount of anti-Growth Hormone polyclonal antibodies, 2) wash away of unbound materials from samples, 3) binding of a second biotinylated anti-Growth Hormone polyclonal antibody to the captured molecules, 4) wash away of unbound materials from samples, 5) conjugation of horseradish peroxidase to the immobilized biotinylated antibodies, 6) wash away of free enzyme conjugates, and 7) quantification of immobilized antibody-enzyme conjugates by monitoring horseradish peroxidase activities in the presence of the substrate 3,3',5,5'-tetramethylbenzidine. The enzyme activity is measured spectrophotometrically by the increased absorbency at 450 nm, corrected from the absorbency at 590 nm, after acidification of formed products. Since the increase in absorbency is directly proportional to the amount of captured rat or mouse Growth Hormone in the unknown sample, the latter can be derived by interpolation from a reference curve generated in the same assay with reference standards of known concentrations of Rat Growth Hormone.

III. REAGENTS SUPPLIED

Each kit is sufficient to run one 96-well plate and contains the following reagents:

A. Microtiter Strip Plate

Coated with affinity purified Polyclonal anti Growth Hormone Antibodies Quantity: 1 plate Preparation: Ready to Use Note: Unused strips should be resealed in the foil pouch with the desiccant provided and stored at 2-8 ℃.

B. Adhesive Plate Sealer

Quantity: 2 sheets Preparation: Ready to Use

III. REAGENTS SUPPLIED (continued)

C. 10X HRP Wash Buffer Concentrate

10X concentrate of 50 mM Tris Buffered Saline containing Tween-20. Quantity: 2 bottles containing 50 mL each Preparation: Dilute 1:10 with distilled or deionized water.

D. Rat/Mouse Growth Hormone Standard

Recombinant Rat Growth Hormone, 0.5 mL/vial, lyophilized. Quantity: 0.5 mL upon hydration Preparation: Reconstitute with 0.5 mL distilled or deionized water.

E. Rat/Mouse Growth Hormone Quality Controls 1 and 2

Rat Growth Hormone, 0.5 mL/vial, lyophilized. Quantity: 0.5 mL/vial upon hydration Preparation: Reconstitute with 0.5 mL distilled or deionized water.

F. Assay Buffer

Buffer containing BSA and 0.08% Sodium Azide Quantity: 40 mL Preparation: Ready to Use

G. Rat/Mouse Growth Hormone Detection Antibody

Pre-titered Biotinylated anti Growth Hormone Polyclonal Antibody Quantity: 12 mL Preparation: Ready to Use

H. Enzyme Solution

Pre-titered Streptavidin-Horseradish Peroxidase Conjugate in Buffer Quantity: 12 mL Preparation: Ready to Use

I. Substrate (Light sensitive, avoid unnecessary exposure to light)

3, 3', 5, 5'-tetramethylbenzidine in buffer Quantity: 12 mL Preparation: Ready to Use.

J. Stop Solution (Caution: Corrosive Solution)

0.3 M HCl Quantity: 12 mL Preparation: Ready to Use

IV. STORAGE AND STABILITY

Prior to use, all components in the kit can be stored up to 2 weeks at 2-8°C. For longer storage (> 2 weeks), freeze Wash Buffer, Assay Buffer, and reconstituted Standards and Controls at $\leq -20^{\circ}$ C. Minimize repeated freeze and thaw of the Growth Hormone Standards and Quality Controls. Refer to expiration dates on all reagents prior to use. Do not mix reagents from different kits unless they have the same lot numbers.

V. REAGENT PRECAUTIONS

A. Sodium Azide

Sodium Azide has been added to certain reagents as a preservative. Although the concentrations are low, Sodium Azide may react with lead and copper plumbing to form highly explosive metal azides. On disposal, flush with a large volume of water to prevent azide build up.

B. Hydrochloric Acid

Hydrochloric Acid is corrosive and can cause eye and skin burns. It is harmful if swallowed and can cause respiratory and digestive tract burns. Avoid contact with skin and eyes. Do not swallow or ingest.

VI. MATERIALS REQUIRED BUT NOT PROVIDED

- 1. Pipettes and Pipette Tips: $10 \ \mu$ L $20 \ \mu$ L or $20 \ \mu$ L $100 \ \mu$ L
- 2. Multi-Channel Pipettes and Pipette Tips: $5 \sim 50 \ \mu L$ and $50 \sim 300 \ \mu L$
- 3. Buffer and Reagent Reservoirs
- 4. Vortex Mixer
- 5. Deionized Water
- 6. Microtiter Plate Reader capable of reading absorbency at 450 nm
- 7. Orbital Microtiter Plate Shaker
- 8. Absorbent Paper or Cloth

VII. SAMPLE COLLECTION AND STORAGE

1. To prepare serum samples, whole blood is directly drawn into a centrifuge tube that contains no anti-coagulant. Let blood clot at room temperature for 30 minutes.

Promptly centrifuge the clotted blood at 2,000 to 3,000 x g for 15 minutes at $4 \pm 2^{\circ}$ C.

Transfer and store serum samples in separate tubes. Date and identify each sample.

Use freshly prepared serum or aliquot and store samples at $\leq -20^{\circ}$ C for later use. For long-term storage, keep at -70°C. Avoid freeze/thaw cycles.

- 2. To prepare plasma samples, whole blood should be collected into centrifuge tubes containing enough K₃EDTA to achieve a final concentration of 1.735 mg/mL and centrifuged immediately after collection. Observe the same precautions in the preparation of serum samples.
- 3. If heparin is to be used as an anticoagulant, the effect on the assay outcome at the dose of heparin used should be pre-determined.
- 4. Avoid using samples with gross hemolysis or lipemia.

VIII. SAMPLE PREPARATION

- 1. No dilution or preparation is needed for normal serum or plasma samples. In the event that any sample is above 50 ng/mL range, dilutions should be performed using the Assay Buffer provided.
- 2. Tissue extracts or cell culture samples may require dilution. Dilutions should be performed using the Assay Buffer provided.

IX. STANDARD AND QUALITY CONTROLS PREPARATION

A. Rat/Mouse Growth Hormone Standard Preparation

- 1. Use care in opening the lyophilized Standard vial. Using a pipette, reconstitute the Rat/Mouse Growth Hormone Standard with 0.5 mL distilled or deionized water into the glass vial to give a 50 ng/mL concentration of Standard. Invert and mix gently, let sit for 5 minutes then mix well.
- 2. Label six tubes 16.7, 5.6, 1.9, 0.62, 0.21, and 0.07 ng/mL. Add 100 μL Assay Buffer to each of the six tubes. Prepare 3 times serial dilutions by adding 50 μL of the 50 ng/mL reconstituted standard to the 16.7 ng/mL tube, mix well and transfer 50 μL of the 16.7 ng/mL standard to the 5.6 ng/mL tube, mix well and transfer 50 μL of the 5.6 ng/mL standard to the 1.9 ng/mL tube, mix well and transfer 50 μL of the 1.9 ng/mL standard to the 0.62 ng/mL tube, mix well and transfer 50 μL of the 0.62 ng/mL standard to the 0.21 ng/mL tube, mix well and transfer 50 μL of the 0.21 ng/mL standard to the 0.07 ng/mL tube, mix well and transfer 50 μL of the 0.21 ng/mL standard to the 0.07 ng/mL tube, mix well and transfer 50 μL of the 0.21 ng/mL standard to the 0.07 ng/mL tube, mix well and transfer 50 μL of the 0.21 ng/mL standard to the 0.07 ng/mL tube, mix well.

Note: Do not use a Repeater pipette. Change tip for every dilution. Wet tip with Standard before dispensing. Unused portions of standard should be stored at \leq -20 °C. Avoid multiple freeze/thaw cycles.

Standard Concentration	Volume of Deionized	Volume of Standard
ng/mL	Water	to Add
	to Add	
50	0.5 mL	0

Standard Concentration ng/mL	Volume of Assay Buffer to Add	Volume of Standard to Add
16.7	100 µL	50 μL of 50 ng/mL
5.6	100 µL	50 μL of 16.7 ng/mL
1.9	100 µL	50 μL of 5.6 ng/mL
0.62	100 µL	50 μL of 1.9 ng/mL
0.21	100 µL	50 μL of 0.62 ng/mL
0.07	100 µL	50 μL of 0.21 ng/mL

IX. STANDARD AND QUALITY CONTROLS PREPARATION (continued)

B. Rat/Mouse Growth Hormone Quality Control 1 and 2 Preparation

1. Use care in opening the lyophilized Quality Control vials. Using a pipette, reconstitute each of the Rat/Mouse Growth Hormone Quality Control 1 and Quality Control 2 with 0.5 mL distilled or deionized water into the glass vials. Invert and mix gently, let sit for 5 minutes then mix well.

X. ASSAY PROCEDURE

Pre-warm all reagents to room temperature prior to setting up the assay.

- 1. Dilute the 10X concentrated Wash Buffer 10 fold by mixing the entire contents of both buffer bottles with 900 mL deionized or distilled water.
- 2. Remove the required number of strips from the Microtiter Assay Plate. Unused strips should be resealed in the foil pouch and stored at 2-8 °C. Assemble the strips in an empty plate holder and wash well with 300 μl of diluted Wash Buffer. Incubate at room temperature for 5 minutes. Decant Wash Buffer and remove the residual amount from all wells by inverting the plate and tapping it smartly onto absorbent towels several times. Do not let wells dry before proceeding to the next step. If an automated machine is used for the assay, use a gentle wash program for all washing steps described in this protocol.
- 3. Add in duplicate, 100 µL Assay Buffer to the blank wells.
- 4. Add in duplicate, 90 μL Assay Buffer to Standard wells, QC1, QC2, and sample wells. (See plate well map for suggested well orientation).
- Add in duplicate, 10 μL Rat/Mouse Growth Hormone Standards in the order of ascending concentration to the appropriate wells. Add in duplicate, 10 μL QC1 and 10 μL QC2 to the appropriate wells. Add sequentially, 10 μL of the unknown samples in duplicate to the remaining wells. For best result all additions should be completed within 30 minutes.
- 6. Cover the plate with plate sealer and incubate at room temperature for 1.5 hours on an orbital microtiter plate shaker set to rotate at moderate speed, approximately 400 to 500 rpm.
- 7. Remove plate sealer and decant solutions from the plate. Tap as before to remove residual solutions in the wells.

X. ASSAY PROCEDURE (continued)

- 8. Wash wells 3 times with diluted Wash Buffer, 300 μ L per well per wash. Decant and tap firmly after each wash to remove residual buffer.
- Add 100 μL Detection Antibody to all wells. Cover the plate with plate sealer and incubate at room temperature for 1 hour on an orbital microtiter plate shaker set to rotate at moderate speed, approximately 400 to 500 rpm.
- 10. Remove plate sealer and decant solutions from the plate. Tap as before to remove residual solutions in the wells.
- 11. Wash wells 3 times with diluted Wash Buffer, 300 μ L per well per wash. Decant and tap firmly after each wash to remove residual buffer.
- 12. Add 100 μ L Enzyme Solution to each well. Cover plate with sealer and incubate with moderate shaking at room temperature for 30 minutes on the microtiter plate shaker.
- 13. Remove sealer, decant solutions from the plate, and tap plate to remove the residual fluid.
- 14. Wash wells 3 times with diluted Wash Buffer, 300 μ L per well per wash. Decant and tap firmly after each wash to remove residual buffer.
- 15. Add 100 μL of Substrate Solution to each well, cover plate with sealer and shake on the plate shaker for approximately 8 to 15 minutes (A longer development time may be needed if using a plate washer). Blue color should be formed in wells of Growth Hormone standards with intensity proportional to increasing concentrations of Growth Hormone.

Note: Please be aware that the color may develop more quickly or more slowly than the recommended incubation time depending on the localized room temperature. Please visually monitor the color development to optimize the incubation time. One can monitor color development using a 370 nm filter, if available, on the spectrophotometer. When the absorbance is between 1.2 and 1.8 at 370 nm, the stop solution can be added to terminate the color development.

16. Remove sealer and add 100 μL Stop Solution [CAUTION: CORROSIVE SOLUTION] and shake plate by hand to ensure complete mixing of solution in all wells. The blue color should turn to yellow after acidification. Read absorbance at 450 nm and 590 nm in a plate reader within 5 minutes and ensure that there are no air bubbles in any well. Record the difference of absorbance units. The absorbance of the highest Growth Hormone standard should be approximately 2.0-3.2, or not to exceed the capability of the plate reader used.

Assay Procedure for Rat/Mouse Growth Hormone ELISA kit (Cat. EZRMGH-45K)

	Step 1	Step 2	Step 3-4	Step 5	Step 6-8	Step 9	Step 9-11	Step 12	Step 12-14	Step 15	Step 15	Step 16	Step 16
Well #	ŗ.		Assay Buffer	Standards/ Controls/ Samples		Detection Ab	Remove residual buffer by ith 300 µl Wash Buffer	Enzyme Solution		Substrate		Stop Solutio	
A1, B1	l Wate	wels	100 µl			100 µl	lual b sh Bu	100 µl		100 µl	re.	100 µl	
C1, D1	Dilute both bottles of 10X Wash Buffer with 900 mL Deionized Water.	h Buffer and incubate for 5 minutes. smartly on absorbent towels	90 µl	10 μl of 0.07 ng/mL Standard	rature.		hour at Room Temperature. Remove residual buffe bsorbent towels. Wash 3X with 300 µl Wash Buffer		at Room Temperature /ash Buffer		Seal, Agitate, Incubate 8-15 minutes at Room Temperature.		
E1, F1	nL Dei	Wash plate with 300 µl Wash Buffer and incubate at room temperature for 5 minutes. residual buffer by tapping smartly on absorbent	90 µl	10 μl of 0.21 ng/mL Standard	ite, Incubate 1.5 hour at Room Temperature. Wash 3X with 300 μl Wash Buffer		Remov th 300		Temp er		n Tem		÷
G1, H1	n 900 n	fer and minut tly on	90 µl	10 μl of 0.62 ng/mL Standard	oom J h Buff		ure. F 3X wi		Room h Buff		t Roon		590 nm.
A2, B2	er with	sh Buf e for 5 J smar	90 µl	10 μl of 1.9 ng/mL Standard	ur at R J Was		nperat Wash		tes at ul Was		utes a		
C2, D2	n Buff	ul Was erature apping	90 µl	10 μl of 5.6 ng/mL Standard	1.5 hoi 1.300 µ		m Ten wels.		minut 300 µ		5 min		uu 05t
E2, F2	K Wasl	h 300 tempe sr by ta	90 µl	10 μl of 16.7 ng/mL Standard	ubate ⁻ 3X with		at Roo ent to		ate 30 3X with		ate 8-1		ice at ²
G2, H2	of 10)	late with 300 μl Wash Buffer and in at room temperature for 5 minutes. Ial buffer by tapping smartly on abs	90 µl	10 μl of 50 ng/mL Standard	Agitate, Incubate Wash 3X wit				, Incub Nash 3		Incuba		Read Absorbance at 450 nm and
A3, B3	oottles	ash pla at esidua	90 µl	10 µl of QC 1	Agita		bate 1 ly on <i>a</i>		gitate		gitate,		ad Abs
C3, D3	both k	Wash plate with 300 µl Was at room temperature Remove residual buffer by tapping	90 µl	10 µl of QC 2	Seal,		e, Incu smart		Seal, Agitate, Incubate 30 minutes at Room Te Wash 3X with 300 µl Wash Buffer		eal, A(Re
E3, F3	Dilute	Ren	90 µl	10 µl of Sample			, Agitate, Incubate 1 tapping smartly on		0,		S		
G3, H3 ↓			90 µl	10 µl of Sample		↓	Seal, , ta	↓		↓ ↓		↓	

XI. MICROTITER PLATE ARRANGEMENT

Rat/Mouse Growth Hormone ELISA

	1	2	3	4	5	6	7	8	9	10	11	12
А	Blank	1.9 ng/mL	QC 1	Sample 3								
в	Blank	1.9 ng/mL	QC 1	Sample 3								
С	0.07 ng/mL	5.6 ng/mL	QC 2	Etc.								
D	0.07 ng/mL	5.6 ng/mL	QC 2									
Е	0.21 ng/mL	16.7 ng/mL	Sample 1									
F	0.21 ng/mL	16.7 ng/mL	Sample 1									
G	0.62 ng/mL	50 ng/mL	Sample 2									
н	0.62 ng/mL	50 ng/mL	Sample 2									

XII. CALCULATIONS

The dose-response curve of this assay fits best to a 5-parameter logistic equation. The results of unknown samples can be calculated with any computer program having a 5-parameter logistic function.


Note: When sample volumes assayed differ from 10 μ L, an appropriate mathematical adjustment must be made to accommodate for the dilution factor (e.g., if 5 μ L of sample is used, then calculated data must be multiplied by 2). When sample volume assayed is less than 10 μ L, compensate the volume deficit with Assay Buffer.

XIII. INTERPRETATION

A. Acceptance Criteria

- 1. The assay will be considered accepted when all Quality Control values fall within the calculated Quality Control Range. If any QC's fall outside the control range, review results with a supervisor.
- 2. If the difference between duplicate results of a sample is >10% CV, repeat the sample.
- 3. The limit of sensitivity of this assay is 0.07 ng/mL Rat/Mouse Growth Hormone (10 μ L sample size).
- 4. The appropriate range of this assay is 0.07 ng/mL to 50 ng/mL Rat/Mouse Growth Hormone (10 μ L sample size). Any result greater than 50 ng/mL in a 10 μ L sample should be diluted using Assay Buffer, and the assay repeated until the results fall within range. Tissue extracts or cell culture media samples greater than 50 ng/mL in a 10 μ L sample should be diluted in Assay Buffer.

XIV. STANDARD CURVE

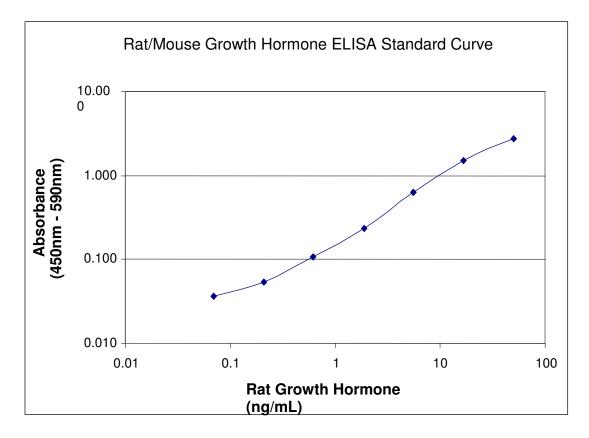

Typical Standard Curve, not to be used to calculate results

XV. ASSAY CHARACTERISTICS

A. Sensitivity

The lowest level of Rat/Mouse Growth Hormone that can be detected by this assay is 0.07 ng/mL when using a 10 μ L sample size.

B. Specificity

Rat Growth Hormone	100%
Rat ACTH	n.d.
Rat TSH	n.d.
Rat FSH	n.d.
Rat LH	n.d.
Rat Prolactin	n.d.
Rat BDNF	n.d.
Human Growth Hormone	n.d.

n.d. = Not Detectable

May cross react with hamster, canine, and primate growth hormone.

C. Precision

Intra-Assay Variation

Sample No.	Mean GH Levels (ng/mL)	Intra-Assay % CV
1	4.1	2.3
2	2.7	4.3
3	2.6	2.1
4	6.1	1.7

Inter-Assay Variation

Sample No.	Mean GH Levels (ng/mL)	Inter-Assay % CV
1	11	4.5
2	5.6	3.2
3	5.2	4.9

The assay variations of Millipore Rat/Mouse Growth Hormone ELISA kits were studied on four rat serum samples with varying concentrations of endogenous Growth Hormone. The mean intra-assay variation was calculated from results of six duplicate determinations in each assay of the indicated samples. The mean inter-assay variation of each sample was calculated from results of four separate assays with duplicate samples in each assay.

XV. ASSAY CHARACTERISTICS (continued)

D. Recovery

Spike & Recovery of Rat Growth Hormone in Rat Serum Samples

Sample No.	GH Added (ng/mL)	Expected (ng/mL)	Observed (ng/mL)	% of Recovery
110.	(119/1112)	(119/1112)	(iig/iiic)	riccovery
1	0	1.74	1.74	100
	3.125	4.865	4.79	98
	6.25	7.99	7.82	98
	12.5	14.24	13.91	98
2	0	3.13	3.13	100
	3.125	6.255	6.13	98
	6.25	9.38	9.15	98
	12.5	15.63	15.78	101
3	0	2.42	2.42	100
	3.125	5.545	5.35	96
	6.25	8.67	8.39	97
	12.5	14.92	14.92	100

Varying amounts of Rat Growth Hormone were added to three rat serum samples and the Growth Hormone content was determined. The % of recovery = observed Growth Hormone concentrations/expected Growth Hormone concentrations x 100%.

XV. ASSAY CHARACTERISTICS (continued)

E. Linearity

Sample	Volume	Expected	Observed	% Of
No.	Sampled	(ng/mL)	(ng/mL)	Expected
	(μL)	((**9,***=)	
1	10	20	20	100
	7.5	15	15.5	103
	5	10	11.2	112
	2.5	5	5.9	118
2	10	4.8	4.8	100
	7.5	3.6	3.6	100
	5	2.4	2.5	104
	2.5	1.2	1.3	108
3	10	3.9	3.9	100
	7.5	2.9	2.8	97
	5	1.95	1.97	101
	2.5	.98	1.0	102

Effect of Serum Dilution

Three rat serum samples with the indicated sample volumes were assayed. Required amounts of matrix were added to compensate for lost volumes below 10 μ L. The resulting dilution factors of 1.0, 1.3, 2.0, and 4.0 representing 10 μ L, 7.5 μ L, 5 μ L, and 2.5 μ L sample volumes assayed, respectively, were applied in the calculation of observed Growth Hormone concentrations. % expected = observed/expected x 100%.

XVI. QUALITY CONTROLS

The ranges for Quality Control 1 and 2 are provided on the card insert or can be located at the Millipore website <u>www.millipore.com</u>.

XVII. TROUBLESHOOTING GUIDE

- 1. To obtain reliable and reproducible results the operator should carefully read this manual and fully understand all aspects of each assay step before attempting to run the assay.
- 2. Throughout the assay the operator should adhere strictly to the procedures with good laboratory practice.
- 3. Have all necessary reagents and equipment ready on hand before starting. Once the assay has been started all steps should be completed with precise timing and without interruption.
- 4. Avoid cross contamination of any reagents or samples to be used in the assay.
- 5. Make sure all reagents and samples are added to the bottom of each well.
- 6. Careful and complete mixing of solutions in the well is critical. Poor assay precision will result from incomplete mixing or cross well contamination due to inappropriate mixing.
- 7. Remove any air bubble formed in the well after acidification of substrate solution because bubbles interfere with spectrophotometric readings.
- 8. Do not let the absorbency reading of the highest standard reach 3.2 units or higher after acidification.
- 9. High absorbance in background or blank wells could be due to 1) cross well contamination by standard solution or sample or 2) inadequate washing of wells with Wash Buffer or 3) overexposure to light after substrate has been added.

XVIII. REPLACEMENT REAGENTS

Reagents	Cat. #
Rat/Mouse Growth Hormone ELISA Plate	EP45
10X HRP Wash Buffer Concentrate	EWB-HRP
Rat/Mouse Growth Hormone Standards	E8045-K
Rat/Mouse Growth Hormone Quality Controls 1 and 2	E6045-K
Assay Buffer	AB-P
Rat/Mouse Growth Hormone Detection Antibody	E1045
Enzyme Solution	EHRP
Substrate	ESS-TMB
Stop Solution	ET-TMB

XIX. ORDERING INFORMATION

A. To place an order:

For USA Customers:

Please provide the following information to our customer service department to expedite your telephone, fax or mail order:

- 1. Your name, telephone and/or fax number
- 2. Customer account number
- 3. Shipping and billing address
- 4. Purchase order number
- 5. Catalog number and description of product
- 6. Quantity and product size

TELEPHONE ORDERS: Toll Free US (866) 441-8400 (636) 441-8400 FAX ORDERS: (636) 441-8050 MAIL ORDERS: Millipore 6 Research Park Drive

St. Charles, Missouri 63304 U.S.A.

For International Customers:

To best serve our international customers, it is Millipore's policy to sell our products through a network of distributors. To place an order or to obtain additional information about Millipore products, please contact your local distributor.

B. Conditions of Sale

All products are for research or manufacturing use only. They are not intended for use in clinical diagnosis or for administration to human or animals. All products are intended for *in vitro* use only.

C. Material Safety Data Sheets (MSDS)

Material safety data sheets for Millipore products may be ordered by fax or phone. See Section A above for details on ordering.