Code No. 27407

Human c-Met Assay Kit - MCM

INTRODUCTION

c-Met, the product of c-met gene has been reported to be the HGF receptor since HGF induces phosphorylation of c-Met by binding specifically to it. It is a heterodimer protein composed of an α -chain that is linked to a β -chain. The β -chain has the intracellular tyrosine kinase domain, the transmembrane domain and the extracellular domain, and it is tied into the α -chain that is the extracellular domain of extracellular domain, and it is the into the d-chain that is the extracellular domain of c-Met. c-Met exists mainly in epidermal cells. It is found in the digestive tract, prostate gland, seminal vesicles, mammary gland, microglia cells, monocytes and macrophages, and more amount in the liver and kidney. It is considered that c-Met tansmits signals resulting from binding to HGF, such as growth, motility and organ formation, in the organs and cells mentioned above. This kit is designed to measure Human c-Met.

| HGE | Product | l inge |
|-----|---------|--------|

| Code No. | Name | Volume |
|----------|--------------------------------------|---------|
| 27401 | Human Activated HGF Assay Kit -MCM | 96 Well |
| 27402 | Human Total HGF Assay Kit - MCM | 96 Well |
| 27403 | Human Activated HGFA Assay Kit - MCM | 96 Well |
| 27404 | Human Total HGFA Assay Kit - MCM | 96 Weil |
| 27405 | Human HAI-1 Assay Kit - MCM | 96 Well |
| 27406 | Human HAI-2 Assay Kit - MCM | 96 Well |
| 27407 | Human c-Met Assay Kit - MCM | 96 Well |

PRINCIPI F

This kit is a solid phase sandwich ELISA using 2 kinds of high specific antibodies. Tetra Methyl Benzidine (TMB) is used as coloring agent (Chromogen). T strength of coloring is in proportion to the quantities of human c-Met.

MEASUREMENT RANGE

0.78 ~ 50 ng/mL

INTENDED USE

- The Human c-Met Assay Kit is a complete kit for the quantitative determination of human c-Met in serum, EDTA-plasma or cell culture media. Determination of human c-Met is affected by the presence of heparin in samples, so please use EDTA-plasma as a sample instead of heparin plasma.

KIT COMPONENT

| 1 | Precoated plate : Ant | i- Human c-Met Mouse IgG MoAb Affinity Purify | 96Well x 1 |
|---|------------------------|-------------------------------------------------|------------|
| 2 | Labeled antibody Co | | |
| | : (30X) HRP conjugated | Anti-Human c-Met Mouse IgG Fab' Affinity Purify | 0.4mL x 1 |
| 3 | Standard | : Human c-Met | 0.5mL x 2 |
| 4 | EIA buffer | | 30mL x 1 |
| 5 | Solution for Labeled | antibody: 1% BSA, 0.05%Tween 20 in PBS | 12mL x 1 |
| 6 | Chromogen | : TMB solution | 15mL x 1 |
| 7 | Stop solution | : 1N H ₂ SO ₄ | 12mL x 1 |
| 8 | Wash buffer Conc. | : (40X) 0.05% Tween20 in phosphate buffer | 50mL x 1 |

OPERATION MANUAL

1. Materials needed but not supplied

| Plate reader (450nm) | · Micropipette and tip |
|--------------------------------------------|-------------------------------------|
| · Graduated cylinder and beaker | Deionized water |
| Incubator (37°C + 1°C) | · Graph paper (log/log |

· Tube for dilution of Standard · Paper towel · Refrigerator (as 4°C) · Washing bottle for precoated plate

· Disposable test tube for "2, Labeled antibody Conc." and "6, Chromogen"

2. Preparation

Preparation of wash buffer 1)

"8, Wash buffer Conc." is a concentrated (40X) buffer. The temperature of "8, Wash buffer Conc." shall be adjusted to room temperature and then, mix it gently and completely before use. Dilute 50mL of *8, Wash buffer Conc." with 1,950mL of deionized water and mix it. This is the wash buffer for use. This prepared wash buffer shall be stored in refrigerator and used within 2 weeks after dilution.

weeks after cliuiton.

Preparation of Labeled antibody

"2, Labeled antibody Conc." is a concentrated (30X). Dilute "2, Labeled antibody Conc." with "5, Solution for Labeled antibody" in 30 times according to required quantity into a disposable test tube. Use this resulting solution as Labeled antibody

Example)

In case you use one slit (8 well), the required quantity of Labeled antibody is $800\,\mu$ L. (Dilute $30\,\mu$ L of "2, Labeled antibody Conc." with $870\,\mu$ L of "5, Solution for Labeled antibody" and mix it. And use the resulting solution by 100 μ L in each well.)

This operation should be done just before the application of Labeled

antibody.
The remaining "2, Labeled antibody Conc." should be stored at 4°C in firmly

Preparation of Standard

Put just <u>0.5 mL</u> of deionized water into the vial of "3, Standard" and mix it gently and completely. This solution is 100 ng/mL Human c-Met standard.

Dilution of Standard


Prepare 8 tubes for dilution of "3, Standard". Put 230 µL each of "4, EIA

Specify the following concentration of each tube

| Tube-1 | 50 ng/mL | |
|--------|------------|--------------------|
| Tube-2 | 25 ng/mL | |
| Tube-3 | 12.5 ng/mL | |
| Tube-4 | 6.25 ng/mL | |
| Tube-5 | 3.13 ng/mL | |
| Tube-6 | 1.56 ng/mL | |
| Tube-7 | 0.78 ng/mL | |
| Tube-8 | 0 ng/mL | (Test Sample Blank |
| | | |

Put 230 $\mu\,\text{L}$ of Standard solution into tube-1 and mix it gently. Then, put 230 $\mu\,\text{L}$ of tube-1 mixture into tube-2. Dilute two times standard solution in series to set up 7 points of diluted standard between 50 ng/mL and 0.78 ng/mL. Tube-8 is the test sample blank as 0 ng/mL.

See following picture

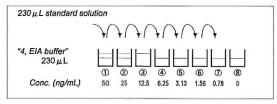

5) Dilution of test sample

Test sample may be diluted with "4, EIA buffer" if the need arises. If the concentration of Human c-Met in samples may not be estimated in advance, the pre-assay with several different dilutions will be recommended to determine the proper dilution of samples.

3. Measurement procedure

All reagents shall be brought to room temperature approximately 30 minutes before use. Then mix it gently and completely before use. Confirm no change in quality of the reagents. Standard curve shall be prepared simultaneously with the measurement of test samples.

| Test Sample | Standard | Test Sample Blank | Reagent Blank |
|------------------------|-----------------------------------------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Test sample 100 μ L | Diluted standard (Tube 1~7) 100 μL | EIA buffer (Tube-8) 100 μL | EIA buffer 100 μL |
| Incubation for 6 | 60 minutes at 37° | C with plate lid | |
| | Washing 7 times | | |
| 100 μ L | 100 μ L | 100 μL | |
| Incubation for | 30 minutes at 4% | C with plate lid | |
| NC | Washing 9 times | 1 | |
| 100 μ L | 100 μ L | 100 μL | 100μL |
| bation for 30 mi | nutes at room te | mperature (shield | led) |
| 100 µ L | 100 µ L | 100 µ L | 100 µ L |
| | Test sample 100 μL Incubation for 6 100 μL Incubation for 100 μL bation for 30 mi | Test sample $100\mu\text{L}$ Diluted standard (Tube 1~7) $100\mu\text{L}$ Incubation for 60 minutes at 37 Washing 7 times $100\mu\text{L}$ $100\mu\text{L}$ Incubation for 30 minutes at 4% Washing 9 times $100\mu\text{L}$ $100\mu\text{L}$ bation for 30 minutes at room te | Test sample 100 μ L Diluted standard (Tube 1~7) 100 μ L EIA buffer (Tube-8) 100 μ L Incubation for 60 minutes at 37°C with plate lid Washing 7 times 100 μ L 100 |

- 1) Determine wells for reagent blank. Put 100 μ L each of "4, EIA buffer" into the
- Determine wells for test sample blank, test sample and diluted standard.
 Then, put 100 μ L each of test sample blank (tube-8), test sample and dilutions of standard (tube-1~7) into the appropriate wells.
 Incubate the precoated plate for 60 minutes at 37°C after covering it with

plate lid.

Wash each well of the precoated plate vigorously with wash buffer using washing bottle. Then, fill each well with wash buffer and leave the precoated washing bottle. Then, fill each well with wash buffer and leave the precoated plate for 15-30 seconds. Remove wash buffer completely from the precoated plate by snapping. This procedure must be repeated more than 7 times. Then, remove the remaining liquid from all wells completely by snapping the Then, remove the fernal liquid from all wells completely by snapping the precoated plate onto paper towel. In case of using plate washer, after 4 times washing with plate washer, washing with above washing bottle must be repeated 3 times. Pipette $100~\mu L$ of labeled antibody solution into the wells of test samples, diluted standard and test sample blank.

Incubate the precoated plate for 30 minutes at 4°C after covering it with plate

Wash the precoated plate 9 times in the same manner above 4).

- *6, Chromogen* should be taken the required quantity into a disposable test tube. Then, pipette 100 μ L from the test tube into the wells. Please avoid to return the rest of test tube into "6, Chromogen" bottle due to avoid to cause of contamination.
- Incubate the precoated plate for 30 minutes at room temperature in the dark.
- The liquid will turn blue by the addition of "6, Chromogen". Pipette $100\,\mu L$ of "7, Stop solution" into the wells. Mix the liquid by tapping the side of precoated plate. The liquid will turn yellow by the addition of "7,
- 11) Remove any dirt or drop of water on the bottom of the precoated plate and confirm there is no bubble on the surface of the liquid. Then, run the plate reader and conduct measurement at 450nm. The measurement shall be done within 30minutes after the addition of "7, Stop solution"



SPECIAL ATTENTION

- Test samples should be measured soon after the collection. In case of the storage of test samples, they should be stored under frozen conditions and do not repeat freeze/thaw cycles. Thaw the test samples at low temperature and mix them completely before measurement.
- Test samples should be diluted with "4, EIA buffer", if the need arises.
- The measurement of test samples and standard in duplicate is
- Use test samples in neutral pH range. The contaminations of organic solvent
- may affect the measurement.

 Use only wash buffer contained in this kit for washing the precoated plate.

 Insufficient washing may lead to the failure in measurement.
- Remove the wash buffer completely by tapping the precoated plate on paper

Do not wipe wells with paper towel.


- "6, Chromogen" should be stored in the dark due to its sensitivity against light.
 "6, Chromogen" should be avoided contact with metals.
- Measurement should be done within 30 minutes after addition of "7, Stop

CALCULATION OF TEST RESULT

Subtract the absorbance of test sample blank from all data, including standards and unknown samples before plotting. Plot the subtracted absorbance of the standards against the standard concentration on log-log graph paper. Draw the best smooth curve through these points to construct the standard curve. Read the concentration for unknown samples from the standard curve.

Example of standard curve

| Conc. (ng/mL) | Absorbance (450nm) |
|--------------------------|-----------------------|
| 50 | 2.549 |
| 25 | 1.199 |
| 12.5 | 0.617 |
| 6.25 | 0.322 |
| 3.13 | 0.158 |
| 1.56 | 0.092 |
| 0.78 | 0.053 |
| 0 (Test Sample Blank) | 0.012 |

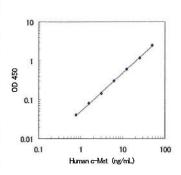

The typical standard curve is shown above. This curve can not be used to derive test results. Please run a standard curve for each assay.

PERFORMANCE CHARACTERISTICS

1. Titer Assay

| Specimen | Titer (X) | Measurement Value (ng/mL) | Theoretical Value (ng/mL) | % |
|----------------------------|--------------|------------------------------|------------------------------|------|
| | 2 | 23.03 | 25.00 | 92.1 |
| 10% FCS added RPMI-1640 | 4 | 10.93 | 12.50 | 87.4 |
| | 8 | 5.74 | 6.25 | 91.8 |
| Human Serum | 2 | 19.81 | 31.82 | 62.3 |
| | 4 | 11.19 | 16.74 | 66.8 |
| | 8 | 6.98 | 9.07 | 77.0 |
| Human Plasma (EDTA) | 2 | 23.83 | 32.50 | 73.3 |
| | 4 | 12.09 | 16.81 | 71.9 |
| | 8 | 7.39 | 9.00 | 82.1 |

2. Added Recovery Assay

| Specimen | Theoretical Value (ng/mL) | Measurement Value (ng/mL) | % |
|---------------------------------|------------------------------|------------------------------|------|
| | 25.09 | 24.41 | 97.3 |
| 10% FCS added RPMI-1640 (x4) | 12.59 | 10.63 | 84.4 |
| 14 Mil 10 10 (X 1) | 6.34 | 5.14 | 81.1 |
| Human Serum (x4) | 15.37 | 9.85 | 64.1 |
| | 9.12 | 6.19 | 67.9 |
| | 5.99 | 4.60 | 76.8 |
| Human Plasma (EDTA) (x4) | 15.45 | 9.85 | 63.8 |
| | 9.20 | 6.56 | 71.3 |
| | 6.08 | 4.82 | 79.3 |

3 Intra - Assay

| Measurement Value (ng/mL) | SD value | CV value (%) | n |
|------------------------------|----------|-----------------|----|
| 17.46 | 0.85 | 4.9 | 24 |
| 6.09 | 0.32 | 5.3 | 24 |
| 2.12 | 0.13 | 6.1 | 24 |

4. Inter - Assay

| Measurement Value (ng/mL) | SD value | CV value (%) | n |
|------------------------------|----------|-----------------|----|
| 18.44 | 1.09 | 5.9 | 36 |
| 6.36 | 0.57 | 9.0 | 36 |
| 2.26 | 0.25 | 11.1 | 36 |

5. Specificity

| Compound | Cross Reactivity | |
|-------------|------------------|--|
| Human c-Met | 100% | |
| Human HGF | ≦0.1% | |
| Human HGFA | ≦0.1% | |
| Human HAI-1 | ≦0.1% | |
| Human HAI-2 | ≦0.1% | |

6. Sensitivity

0.26 ng/mL

0.26 righth.

The sensitivity for this kit was determined using the guidelines under the National Committee for Clinical Leboratory Standards (NCCLS) Evaluation Protocols. (National Committee for Clinical Laboratory Standards Evaluation Protocols, SC1, (1989) Villanova, PA: NCCLS.)

- PRECAUTION FOR INTENDED USE AND/OR HANDLING

 1. All reagents should be stored at 2~8°C. All reagents shall be brought to room temperature approximately 30 minutes before use.

 2. "3, Standard" is lyophilized products. Be careful to open this vial.

 3. "7, Stop solution" is a strong acid substance. Therefore, be careful not to contact your skin and clothes with "7, Stop solution" and pay attention to the disposal of "7, Stop solution".

 4. "1, Precoated plate" and "3, Standard" contain sodium azide. Therefore, dispose these materials after diluting them with large quantity of water to avoid the production of explosive metallic azide.
- the production of explosive metallic azide.
- The precipitation may grow in "2, Labeled antibody Conc.", however, there is no problem in the performance.

 Wash hands after handling reagents.

 Do not mix the reagents with the reagents from different lot or different kit.

Do not use the reagents expired.

This kit is for research purpose only. Do not use for clinical diagnosis.

STORAGE AND THE TERM OF VALIDITY

Storage Condition

:2~8℃

The term of validity : 12 months

(The expiry date is specified in outer box.)

REFERENCE

- Kataoka H, Hamasuna R, Itoh H, Kitamura N, Koono M. Activation of hepatocyte growth factor/scatter factor in colorectal carcinoma. Cancer Res. 2000 Nov 1;60(21):6148-59.
- Shimomura T, Denda K, Kitamura A, Kawaguchi T, Kito M, Kondo J, Kagaya S, Qin L, Takata H, Miyazawa K, Kitamura N. Hepatocyte growth factor activator inhibitor, a novel Kunitz-type serine protease inhibitor. J Biol Chem. 1997 Mar 7:272(10):6370-6.
- Kawaguchi T, Qin L, Shimomura T, Kondo J, Matsumoto K, Denda K, Kitamura N. Purification and cloning of hepatocyte growth factor activator inhibitor type 2, a Kunitz-type serine protease inhibitor. J Biol Chem. 1997 Oct 31;272(44):27558-64.

Version 061204 Established Changed product name

<Distributed by> SCETI K.K. DF Kasumigaseki Place, 3-6-7, Kasumigaseki, Chiyoda-ku Tokyo 100-0013 Japan TEL: +81-(0)3-5510-2347

FAX: +81-(0)3-5510-0134 URL: http://www.sceti.co.jp/export/

e-mail: exp-pet@sceti.co.jp

