Code No. 27776 Human APP βCTF Assay Kit - IBL

INTRODUCTION

Alzheimer's disease (AD) was first reported by A. Alzheimer, a German neuropathologist in 1907 and is considered as a major factor of senile dementia. It is known that Amyloid β (A β ; which is a major constituent of senile plaque) is cleaved from Amyloid Precursor Protein (APP; which exists in three main isoforms, APP695, APP751, and APP770) by β -secretase and subsequent γ -secretase (ref. 1). On the one hand, it has been also reported there is β CTF (β -secretase C-terminal fragment), fragments which are cleaved by β-secretase but not by v-secretase

It is thought that measuring β CTF, in addition to A β , soluble APP α and APP β , also

has significance in research of the metabolic pathway of APP. This ELISA can measure human APP βCTF in samples such as cultured cell lysate .

PRINCIPLE

This kit is a solid phase sandwich ELISA using 2 kinds of high specific antibodies. Tetra Methyl Benzidine (TMB) is used as a coloring agent (Chromogen). The strength of coloring is proportional to the quantities of Human APP β CTF.

MEASUREMENT RANGE

0.19 - 12 pmol/L

INTENDED USE

This IBL's assay kit is capable for the measurement of Human APP BCTF in cell lvsate

- Preparation of the cultured cell lysate
- Add IBLysis-I (IBL Lysate buffer #19022) to the cells pellet (2x10⁷ indication).
 Mix thoroughly with a mixer like Vortex.
- 3 Solubilize by rotation at 2-8°C for 30 minutes
- Centrifuge at 10,000 rpm for 10 minutes at 2-8°C.
 Apply the supernatant diluted by "4, EIA buffer" as necessary.

KIT COMPONENT

1	Precoated plate : A	nti-APP-C Rabbit IgG Affinity Purify	96Well x 1
2	Labeled antibody Co	onc.	
	: (30X) HRP conjugated	Anti- Human Aβ (N) (82E1) Mouse IgG MoAb Affinity Purify	0.4mL x 1
3	Standard : Synthet	ic peptide corresponding to Human APP βCTF	0.5mL x 2
4	EIA buffer : 1% BS/	A, 0.05% Tween20 in PBS	30mL x 1
5	Solution for Labeled	antibody : 1% BSA, 0.05% Tween20 in PBS	12mL x 1
6	Chromogen	: TMB solution	15mL x 1
7	Stop solution	: 1N H ₂ SO ₄	12mL x 1
8	Wash buffer Conc.	: (40X) 0.05% Tween20 in phosphate buffer	50mL x 1
OPE	ERATION MANUAL		

Interials needed but not supplied

waterials	needed	ραι ποι	supplied
 Plat 	e reader	(450nm	1)

Refrigerator (as 4°C)

 Plate reader (450nm) 	 Micropipette and tip
 Graduated cylinder and beaker 	 Deionized water

- · Deionized water
 - · Graph paper (log/log)
- Paper towel
- Tube for dilution of Standard · Washing bottle for precoated plate
- · Disposable test tube for "2, Labeled antibody Conc." and "6, Chromogen"

2. Preparation

1)

Preparation of wash buffer "8, Wash buffer Conc." is a concentrated (40X) buffer. Adjust the temperature of "8, Washing buffer Conc." to room temperature and then, mix it gently and completely before use. Dilute 50 mL of "8, Wash buffer Conc." with 1,950 mL of deionized water and mix it. This is the wash buffer for use. This prepared wash buffer shall be stored in refrigerator and used within 2 weeks after dilution.

Preparation of Labeled antibody 2)

"2, Labeled antibody Conc." is a concentrated (30X). Dilute "2, Labeled antibody Conc." with "5, Solution for Labeled antibody" in 30 times according to required quantity into a disposable test tube. Use this resulting solution as Labeled antibody. Example)

In case you use one strip (8 well), the required quantity of Labeled antibody is 800 μ L. (Dilute 30 μ L of "2, Labeled antibody Conc." with 870 μ L of "5, Solution for Labeled antibody" and mix it. And use the resulting solution by 100 µL in each well.)

This operation should be done just before the application of Labeled antibody.

The remaining "2, Labeled antibody Conc." should be stored at 4°C in firmly sealed vial.

Preparation of Standard 3)

Put just 0.5 mL of deionized water into the vial of "3, Standard" and mix it gently and completely. This solution is 24 pmol/L human APP β CTF standard.


Dilution of Standard Prepare 8 tubes for dilution of "3, Standard". Put 230 µL each of "4, EIA buffer" into the tube.

Specify the following concentration of each tube.'

Tube-1	12 pmol/L	
Tube-2	6 pmol/L	
Tube-3	3 pmol/L	
Tube-4	1.5 pmol/L	
Tube-5	0.75 pmol/L	
Tube-6	0.38 pmol/L	
Tube-7	0.19 pmol/L	
Tube-8	0 pmol/L	(Test Sample Blank)

Put 230 µL of Standard solution into tube-1 and mix it gently. Then, put 230 µL of tube-1 mixture into tube-2. Dilute two times standard solution in series to set up 7 points of diluted standard between 12 pmol/L and 0.19 pmol/L. Tube-8 is the test sample blank as 0 pmol/L.

See following picture.

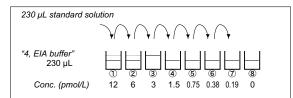

5) Dilution of test sample

Test sample should be diluted with "4, EIA buffer" as necessary. If the quantity of human APP β CTF in samples may not be estimated in advance, the pre-assay with several different dilutions will be recommended to determine the proper dilution of samples

3. Measurement procedure

All reagents shall be brought to room temperature approximately 30 minutes before use. Then mix it gently and completely before use. Make sure of no change in quality of the reagents. Standard of simultaneously with the measurement of test samples. Standard curve shall be prepared

	Test Sample	Standard	Test Sample Blank	Reagent Blank
Reagents	Test sample 100 µL	Diluted standard (Tube 1~7) 100 µL	EIA buffer (Tube-8) 100 µL	EIA buffer 100 μL
	Incubation	overnight at 4°C	with plate lid	
		Washing 7 times	5	
Labeled Antibody 100 µL 100 µL 100 µL				
	Incubation for	60 minutes at 4°	C with plate lid	
		Washing 9 times	5	
Chromogen	100 µL	100 µL	100 µL	100 µL
Incubation for 30 minutes at room temperature (shielded)				
Stop solution	100 µL	100 µL	100 µL	100 µL
Read the plate at 450nm against a Reagent Blank within 30 minutes after addition of Stop solution.				

- 1) Determine wells for reagent blank. Put 100 µL each of "4, EIA buffer" into the
- Determine wells for test sample blank, test sample and diluted standard. 2)
- Determine wells for test sample blank, test sample and diluted standard. Then, put 100 μ L each of test sample blank (tube-8), test sample and dilutions of standard (tube-17) into the appropriate wells. Incubate the precoated plate overnight at 4°C after covering it with plate lid. Wash each well of the precoated plate vigorously with wash buffer using the washing bottle. Then, fill each well with wash buffer and leave the precoated plate laid for 15-30 seconds. Remove wash buffer completely from the precoated plate by snapping. This procedure must be repeated more than 7 times. Then, remove the remaining liquid from all wells completely by snapping the precoated plate onto paper towel. In case of using a plate washer, after 4 times washing with plate washer, washing with above washing bottle must be repeated 3 times. Pipette 100 μ L of labeled antibody solution into the wells of test samples,
- 5)
- Pipette 100 μ L of labeled antibody solution into the wells of test samples, diluted standard and test sample blank.
- 6) Incubate the precoated plate for 60 minutes at 4°C after covering it with plate lid.
- IId. Wash the precoated plate 9 times in the same manner as 4). Take the required quantity of "6, Chromogen" into a disposable test tube. Then, pipette 100 μ L from the test tube into the wells. Please do not return the rest of the test tube to "6, Chromogen" bottle to avoid contamination. Incubate the precoated plate for 30 minutes at room temperature in the dark. The liquid will turn blue by addition of "6, Chromogen". Pipette 100 μ L of "7, Stop solution" into the wells. Mix the liquid by tapping the side of precoated plate. The liquid will turn yellow by addition of "7, Stop solution" 8)
- 9)
- 10) solution'
- Remove any dirt or drop of water on the bottom of the precoated plate and confirm there is no bubble on the surface of the liquid. Then, run the plate 11) reader and conduct measurement at 450 nm against a reagent blank. The measurement shall be done within 30 minutes after addition of "7, Stop solution".

SPECIAL ATTENTION

- 1) Test samples should be measured soon after collection. For the storage of test samples, store them frozen and do not repeat freeze/thaw cycles. Thaw the test samples at a low temperature and mix them completely before measurement.
- Test samples should be diluted with "4, EIA buffer", if the need arises
- 3)
- Duplicate measurement of test samples and standard is recommended. Use test samples in neutral pH range. The contaminations of organic solvent 4) may affect the measurement. Use only wash buffer contained in this kit for washing the precoated plate.
- 5) Insufficient washing may lead to the failure in measurement.
- Remove the wash buffer completely by tapping the precoated plate on paper towel. Do not wipe wells with paper towel. 6)
- 7) "6, Chromogen" should be stored in the dark due to its sensitivity against light.




Instructions Code No. 27776

"6. Chromogen" should be avoided contact with metals Measurement should be done within 30 minutes after addition of "7, Stop solution"

CALCULATION OF TEST RESULT

Subtract the absorbance of test sample blank from all data, including standards and unknown samples before plotting. Plot the subtracted absorbance of the standards against the standard concentration on log-log graph paper. Draw the best smooth curve through these points to construct the standard curve. Read the concentration for unknown samples from the standard curve.

Example of standard curve

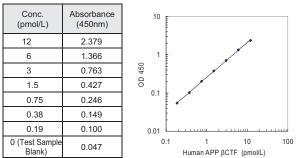

The typical standard curve is shown above. This curve can not be used to * derive test results. Please run a standard curve for each assay.

PERFORMANCE CHARACTERISTICS

1. Titer Assay (Samples with standard added are used.)

Specimen	Titer (X)	Measurement Value (pmol/L)	Theoretical Value (pmol/L)	%
	2	3.25	3.77	86.2
NB-1 Cell lysate	4	1.63	1.94	84.0
(IBLysis-I)	8	0.89	0.99	89.9
	16	0.51	0.47	108.5
Neuro2a	4	1.23	1.58	77.8
Cell lysate	8	0.70	0.80	87.5
(IBLysis-I)	16	0.36	0.41	87.8

IBLysis-I: Lysate buffer by IBL (# 19022)

2. Added Recovery Assay

Specimen	Theoretical Value (pmol/L)	Measurement Value (pmol/L)	%
	3.00	2.92	97.3
IBLysis-I (x2)	1.50	1.41	94.0
	0.75	0.71	94.7

3. Intra - Assay

Measurement Value (pmol/L)	SD value	CV value (%)	n
5.68	0.11	1.9	24
1.38	0.06	4.3	24
0.42	0.02	4.8	24

4. Inter - Assay

[Measurement Value (pmol/L)	SD value	CV value (%)	n
	5.39	0.23	4.3	5
	1.32	0.05	3.8	5
Ī	0.44	0.01	2.3	5


5. Specificity

Compound	Cross Reactivity
Human APP βCTF	100%
Human sAPPβ-Wild Type	≦0.1%
Human sAPPβ-Swedish Type	≦0.1%
Human sAPPα	≦0.1%

6. Sensitivity

0.02 pmol/L The sensitivity for this kit was determined using the guidelines under the National Committee for Clinical Laboratory Standards (NCCLS) Evaluation Protocols. (National Committee for Clinical Laboratory Standards Evaluation Protocols, SC1, (1989) Villanova, PA: NCCLS.)

7. Example of an experiment

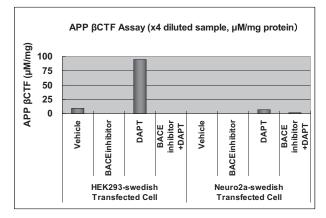

Data provided by Dr. Nobumasa Takasugi, Department of Neuropathology and Neuroscience, Graduate School of Pharmaceutical Sciences, University of Tokyo, Tokyo, Japan

PRECAUTION FOR INTENDED USE AND/OR HANDLING

- All reagents should be stored at 2 8°C. All reagents shall be brought to room 1. temperature approximately 30 minutes before use
- 2
- "3, Standard" is lyophilized products. Be careful to open this vial. "7, Stop solution" is a strong acid substance. Therefore, be careful not to have your skin and clothes contact "7, Stop solution" and pay attention to the disposal of "7, Stop solution". 3.
- Dispose used materials after rinsing them with large quantity of water. Precipitation may occur in "2, Labeled antibody Conc.", however, there is no 5.
- problem in the performance. Wash hands after handling reagents. 6.
- 7 Do not mix the reagents with the reagents from a different lot or kit. Do not use expired reagents.
- 8.
- 9 This kit is for research purpose only. Do not use for clinical diagnosis.

STORAGE AND THE TERM OF VALIDITY

Storage Condition : 2 - 8°C The expiry date is specified on outer box.

REFERENCE

Selkoe DJ. Normal and abnormal biology of the beta-amyloid precursor protein. Annu Rev Neurosci. 1994;17:489-517.

Version 1.3

Amyloid β and APP related products

Code No.	Product Name	Volume	
27711	Human Amyloidβ (1-42) Assay Kit - IBL	96 Well	
27712	Human Amyloidβ (1-42) (N) Assay Kit - IBL	96 Well	
27713	Human Amyloidβ (1-40) Assay Kit - IBL	96 Well	
27714	Human Amyloidβ (1-40) (N) Assay Kit - IBL	96 Well	
27718	Human Amyloidβ (1-40) (FL) Assay Kit - IBL	96 Well	
27720	Mouse/Rat Amyloidβ (1-40) High Specific Assay Kit - IBL	96 Well	
27716	Human Amyloidβ (N3pE-42) Assay Kit - IBL	96 Well	
27418	Human Amyloidβ (N3pE-40) Assay Kit - IBL	96 Well	
27729	Human Amyloidβ (1- x) Assay Kit - IBL	96 Well	
27724	Human sAPPα Assay Kit - IBL	96 Well	
27415	Mouse/Rat sAPPα Assay Kit - IBL	96 Well	
27722	Human sAPPβ-Wild Type Assay Kit - IBL	96 Well	
27723	Human sAPPβ-Swedish Type Assay Kit - IBL	96 Well	
27776	Human APP βCTF Assay Kit - IBL	96 Well	
19022	IBLysis-I (Lysate buffer)	50 mL	

Made in Japan.

<Distributed by>



Tel: +81-3-5510-2347

SCETI K.K. 3-6-7, Kasumigaseki, Chiyoda-ku, Tokyo, 100-0013 Japan

URL: www.sceti.co.jp/export/ e-mail: exp-pet@sceti.co.jp Fax: +81-3-5510-0133